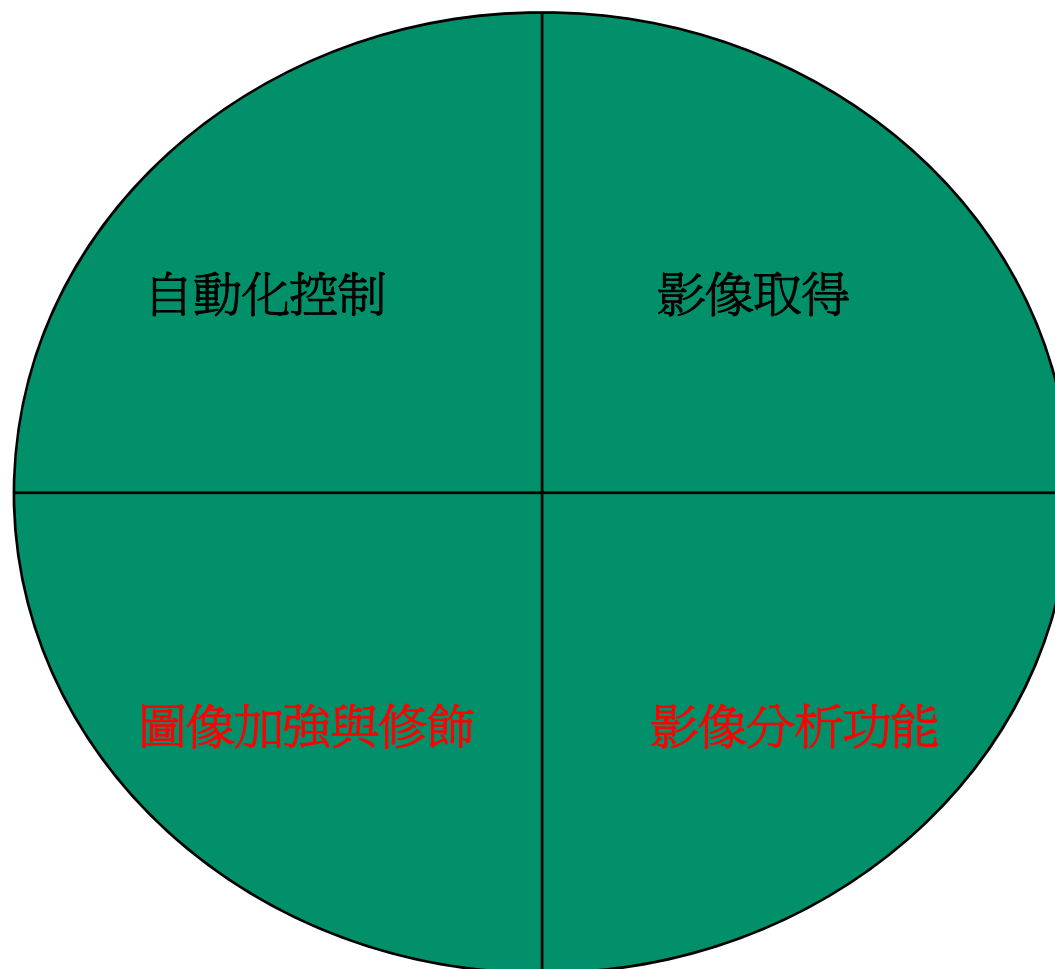
MetaMorph – The Most Powerful Imaging Tool

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
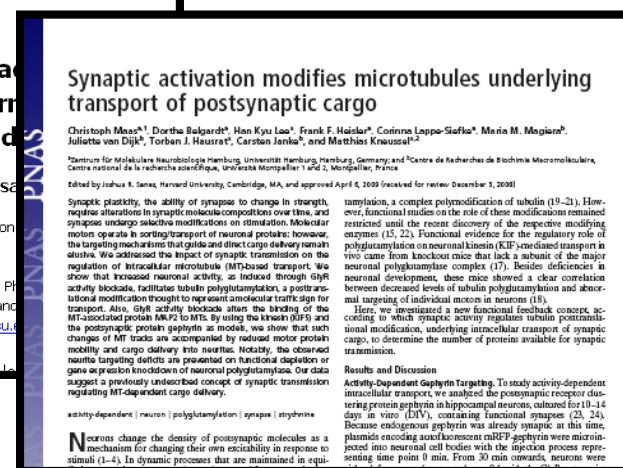
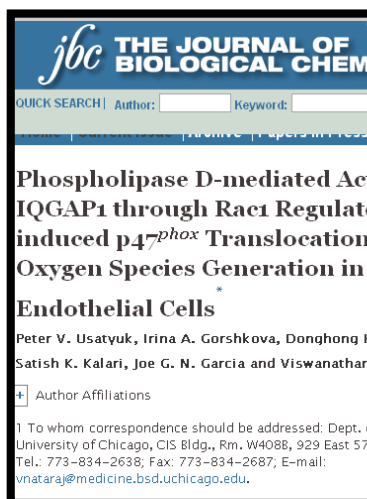
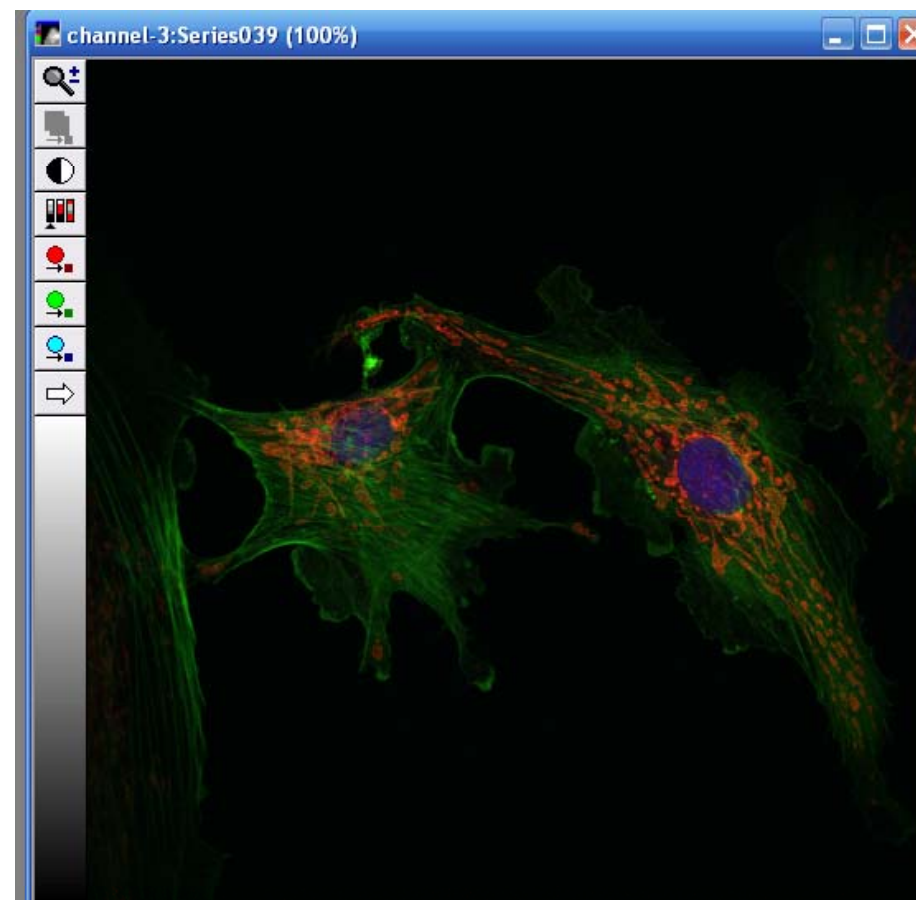
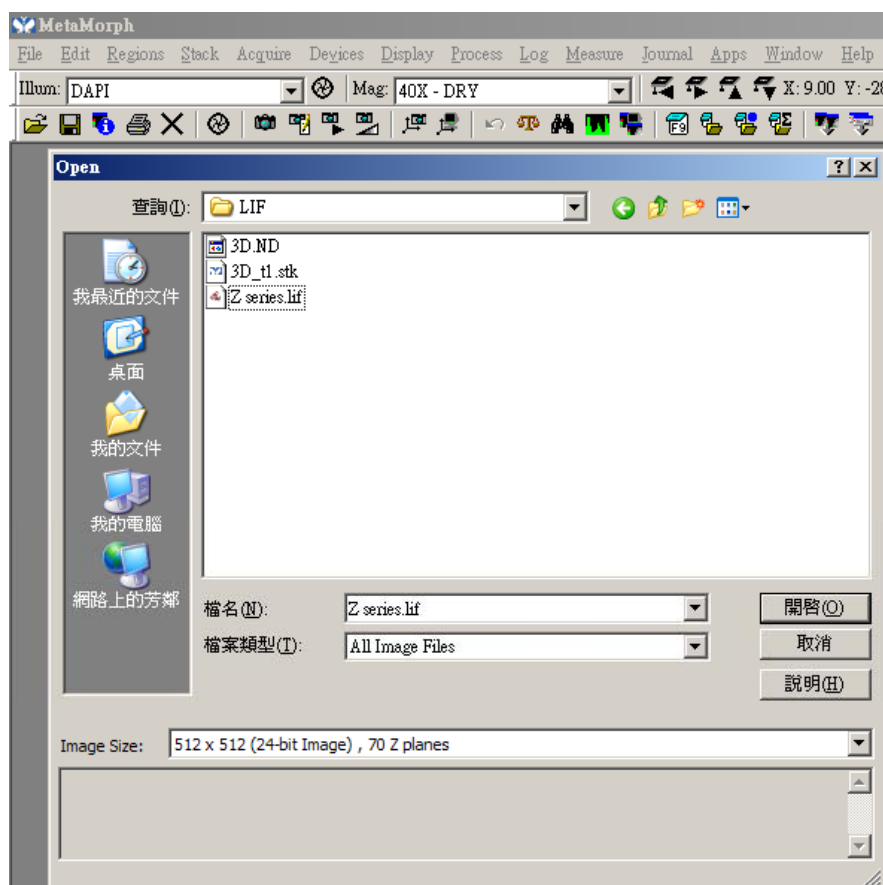
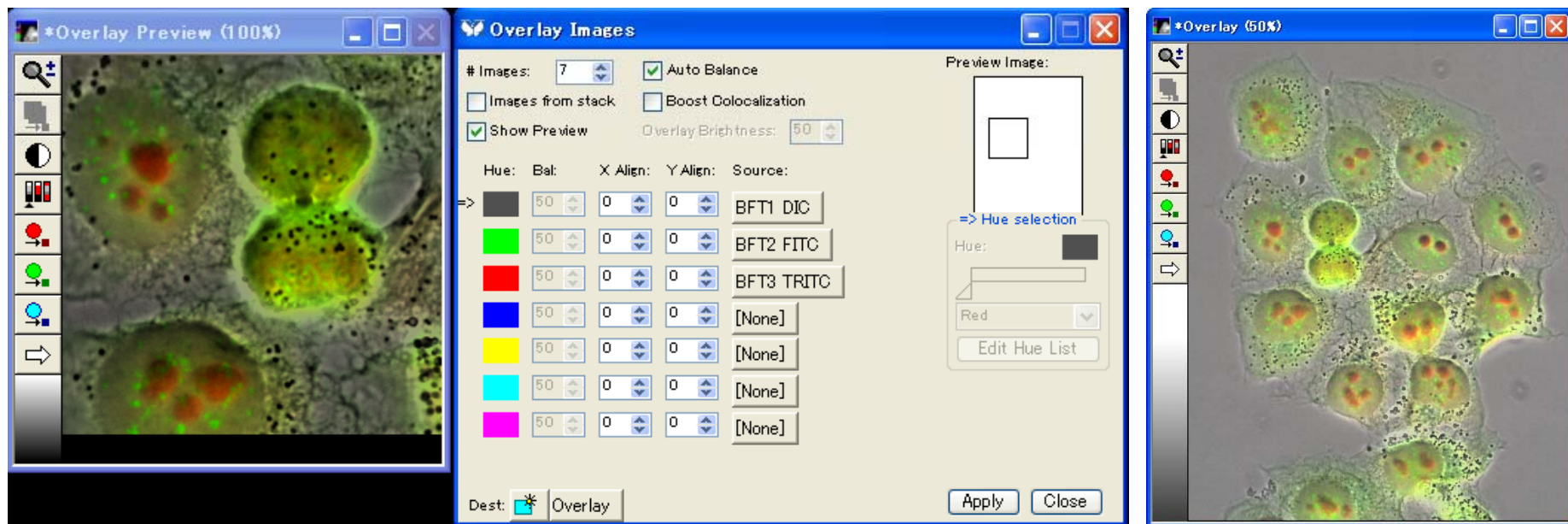
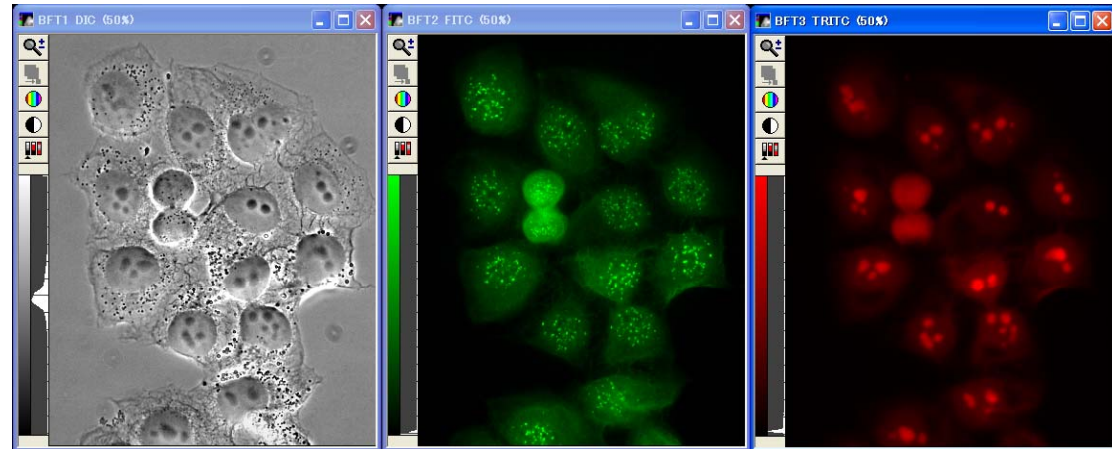
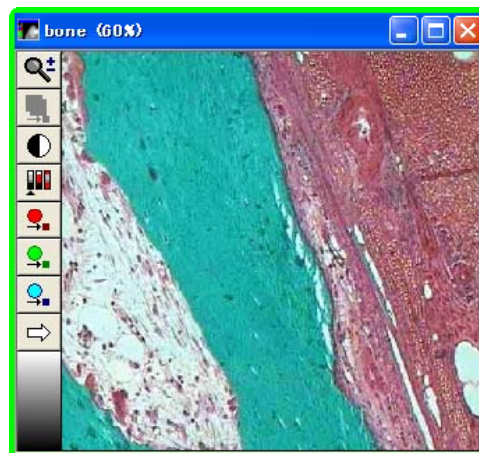
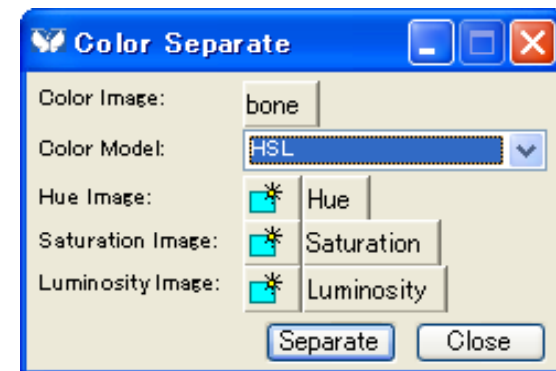
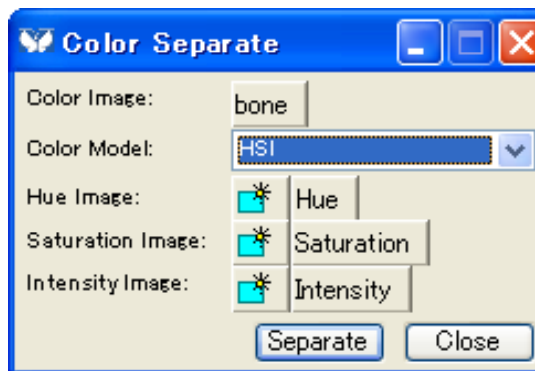
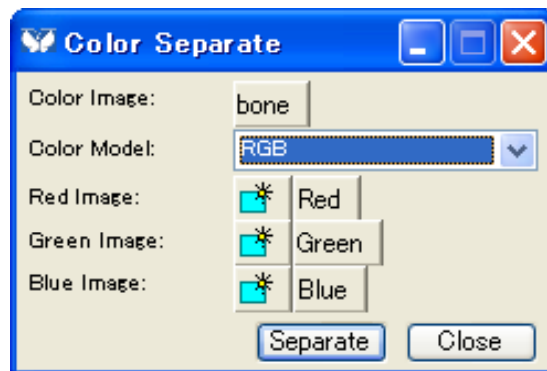
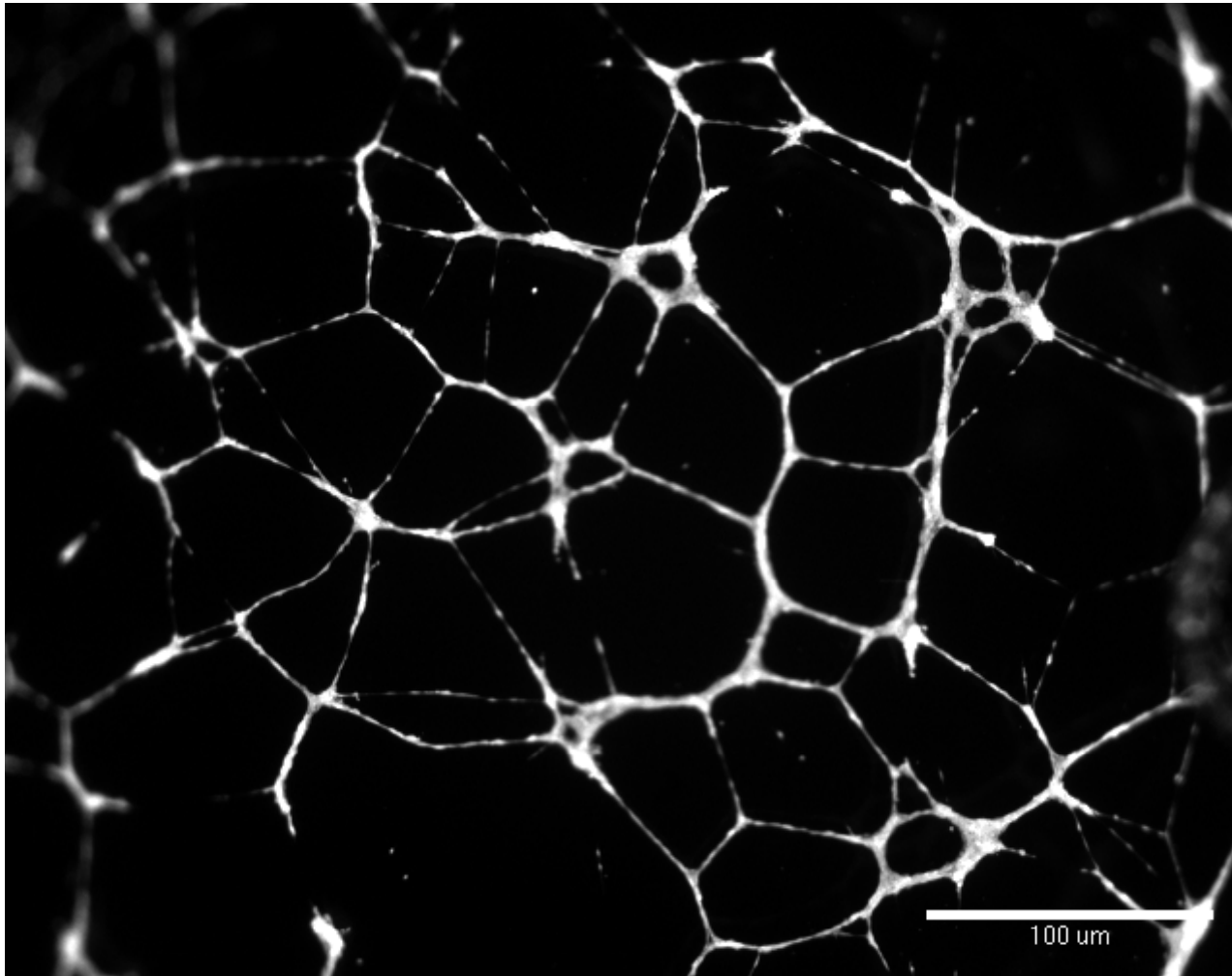
Basic Measurements



Region Measurements (手動量測)

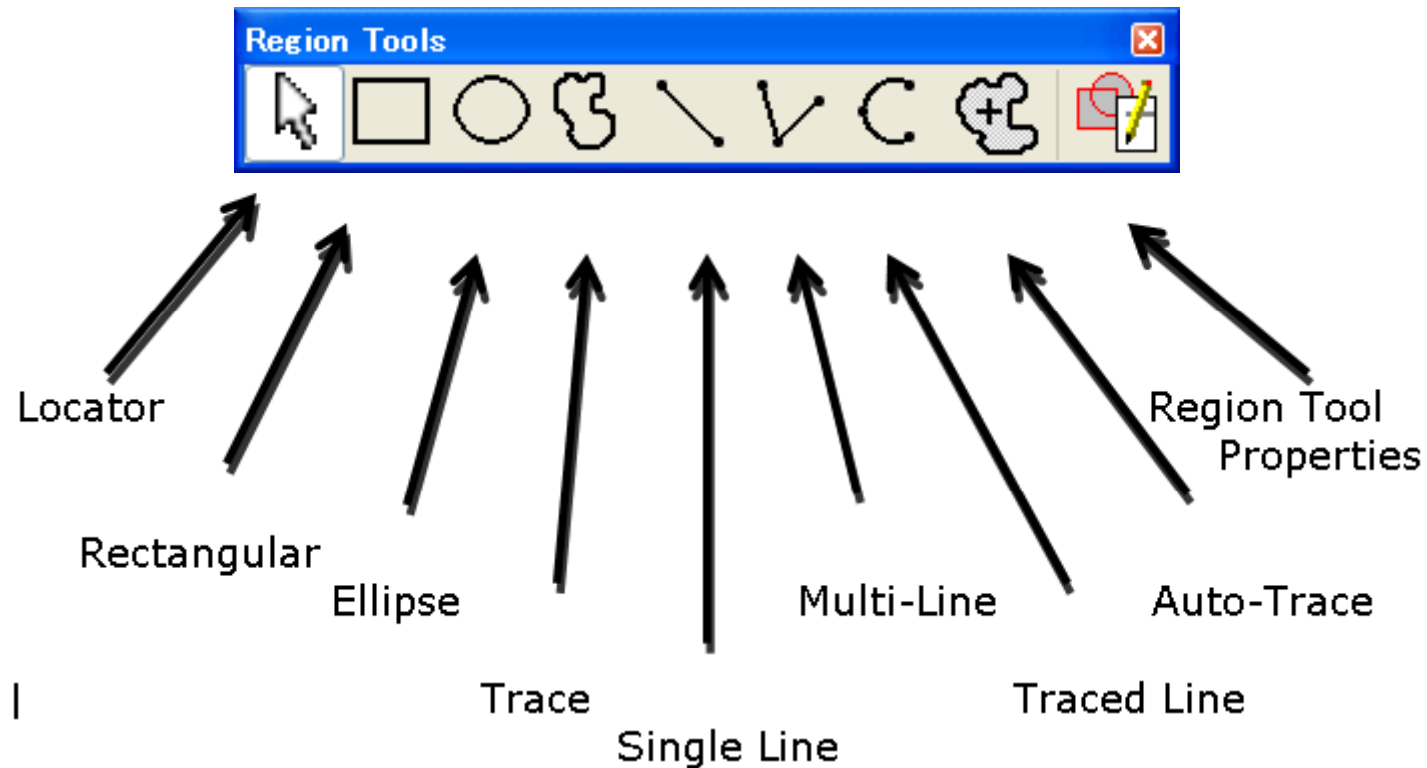➤ Area, Distance, Angle, X/Y Coordinate, Width, Height, Intensity

Region Tooltips Info

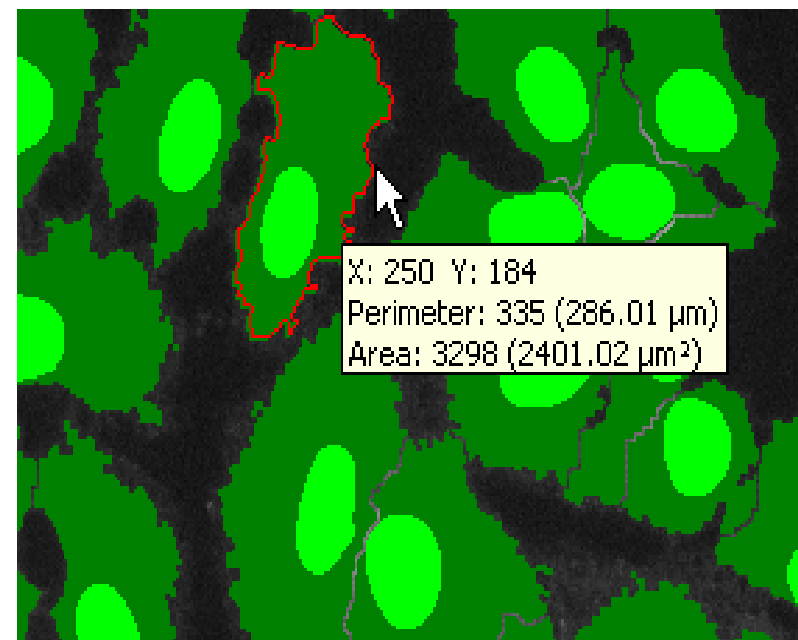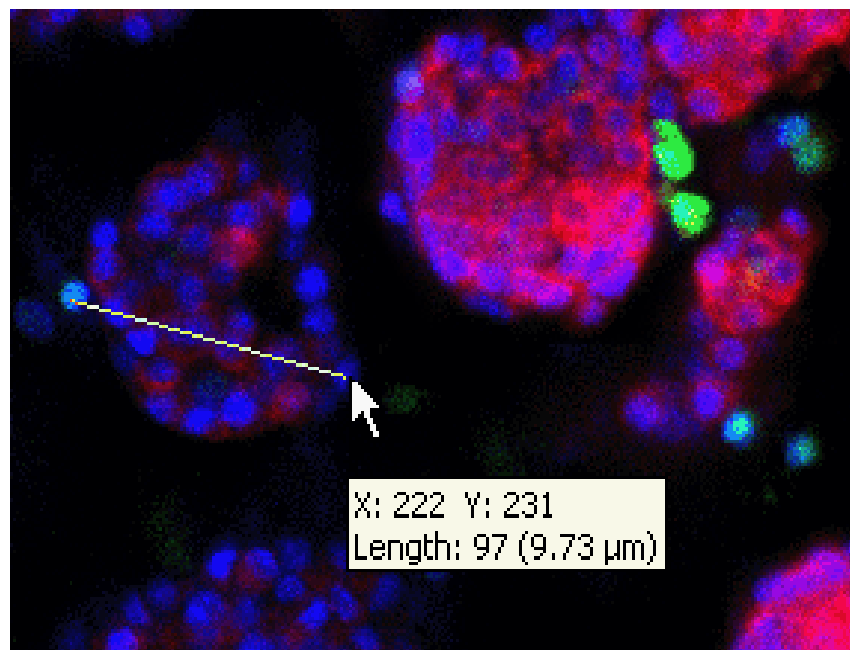Region Measurements - Results


Region Measurements

timelp  Open Log


Include: All Regions  Close


Measurements | Graph | Configure | Labels

Region Label	Image Plane	Average Intensity	Minimum Intensity	Maximum Intensity
1	1	89.8197	8	255
1	2	91.082	12	186
1	3	90.377	10	173
1	4	91.4918	0	185
1	5	93.0656	2	191
1	6	93.6066	4	187
1	7	93.9344	5	185
1	8	95.3115	2	201
1	9	94.5246	8	177
1	10	91.3443	29	145
1	11	89.1803	57	147
1	12	88.4918	59	148
Integrated	-	1989.93	752	3648

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
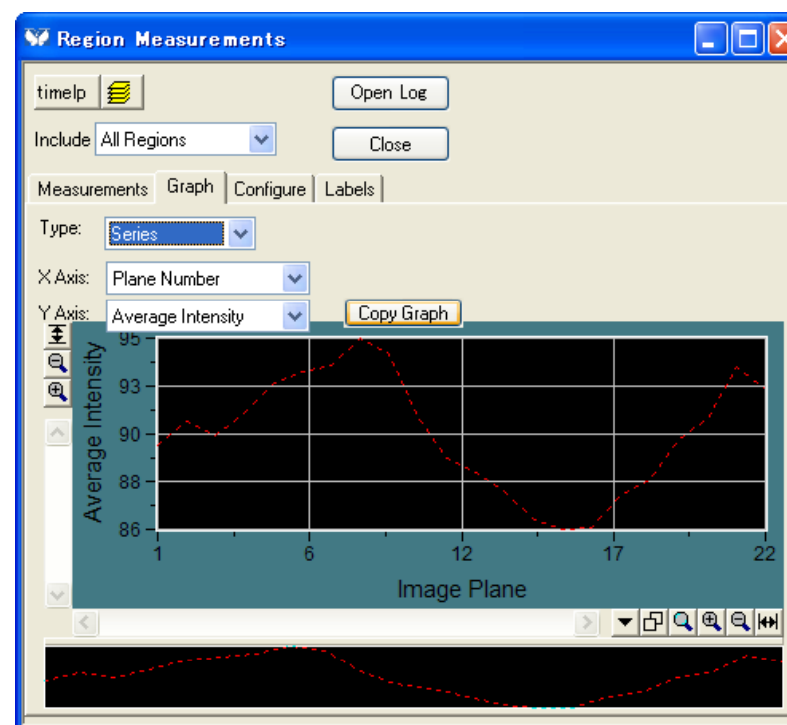
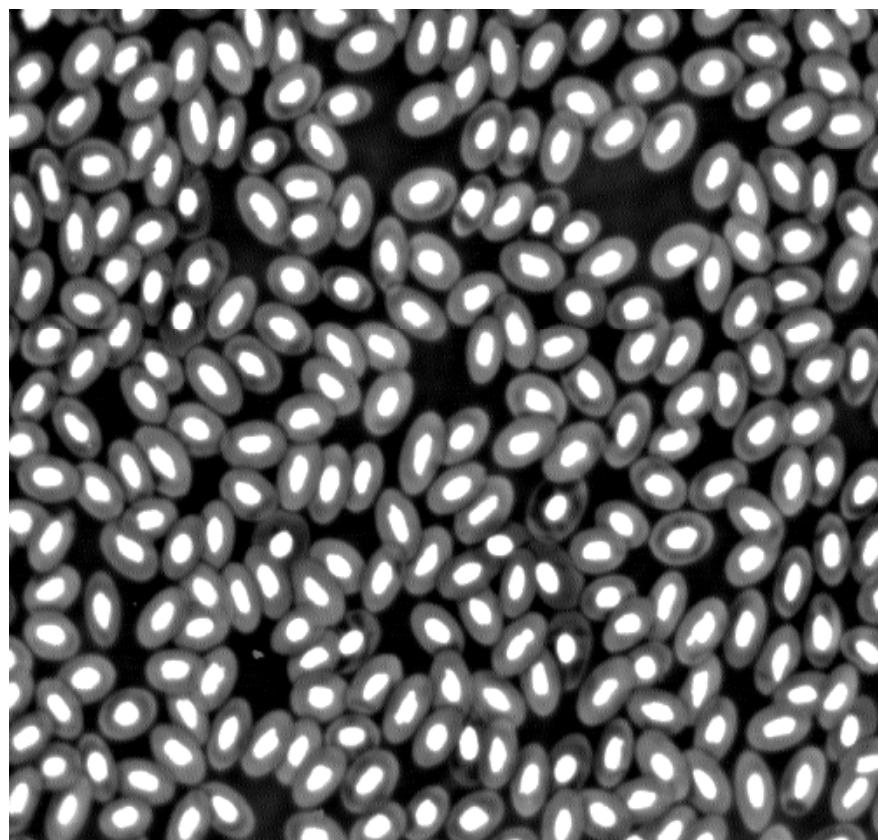
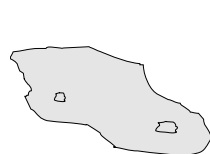
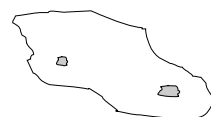
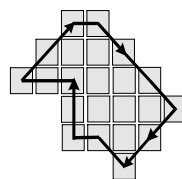
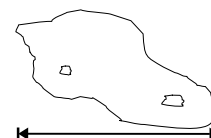
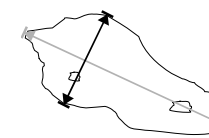
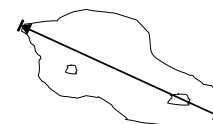
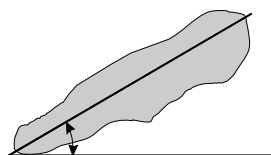
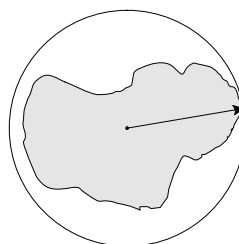
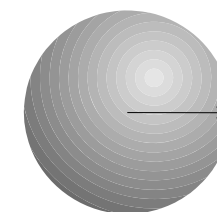
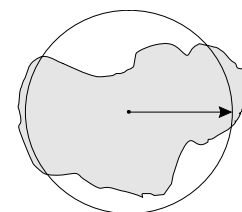
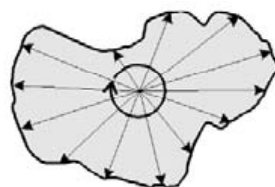
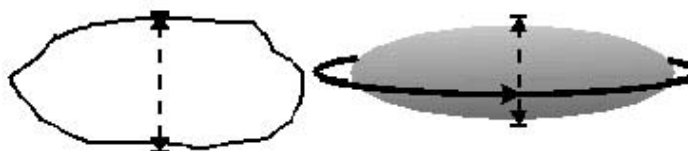
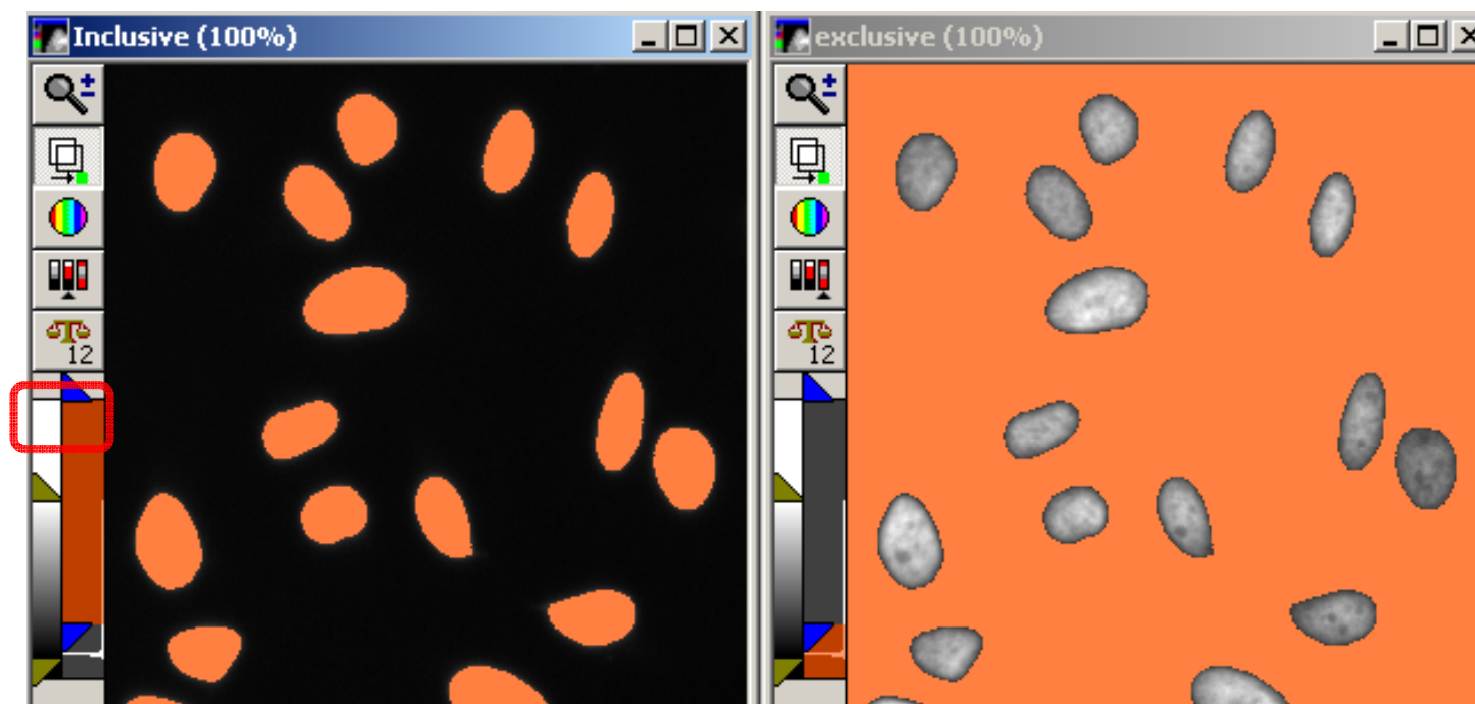
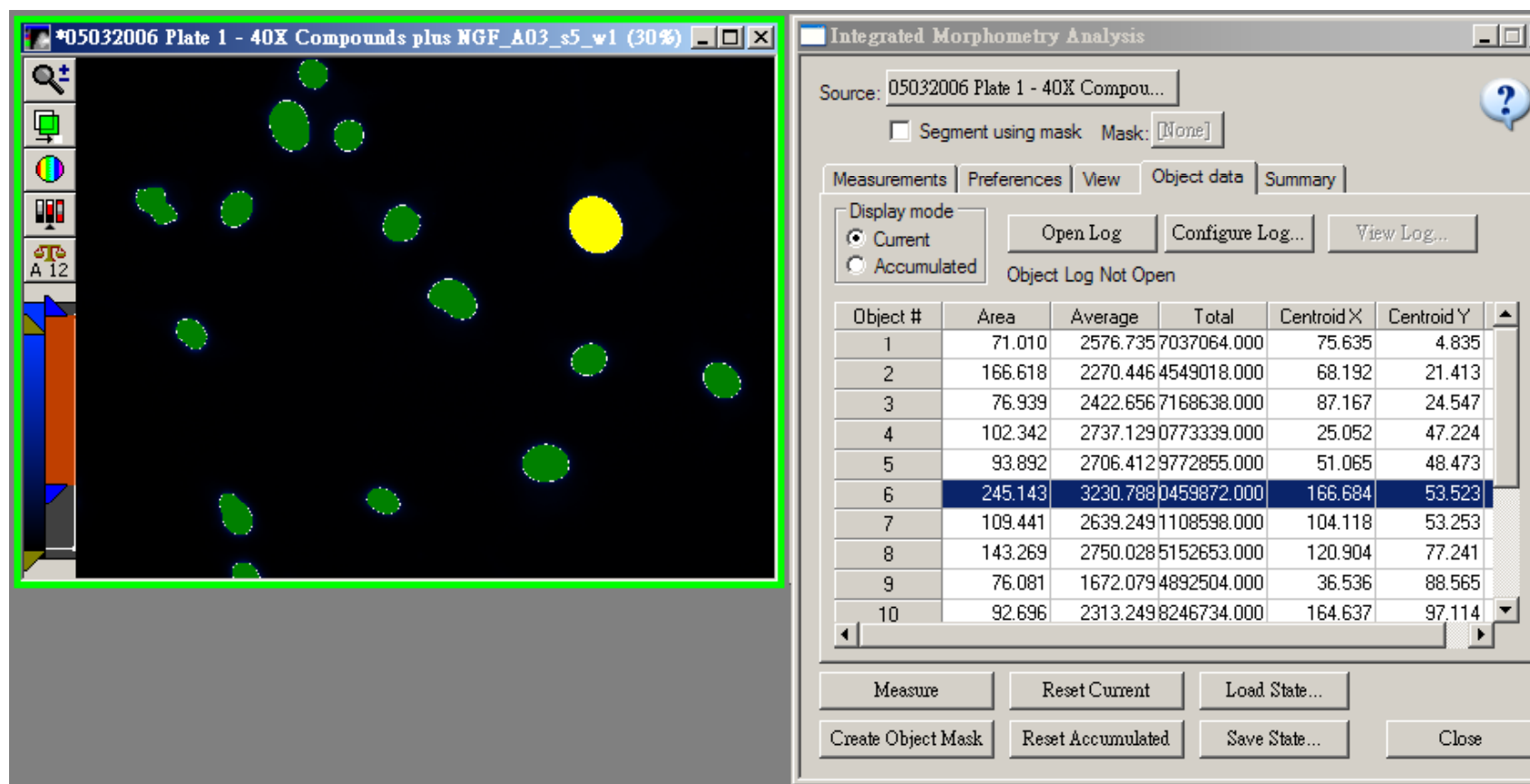
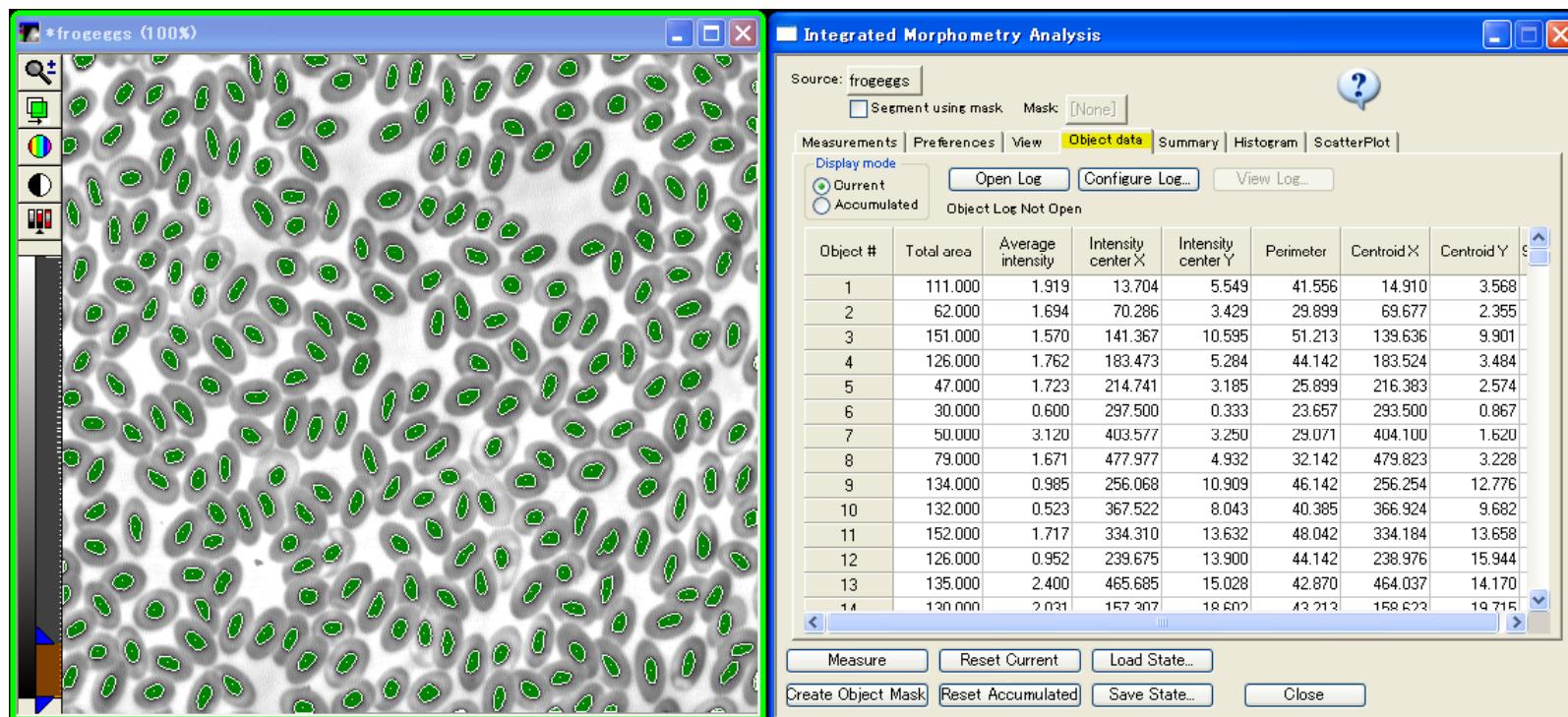
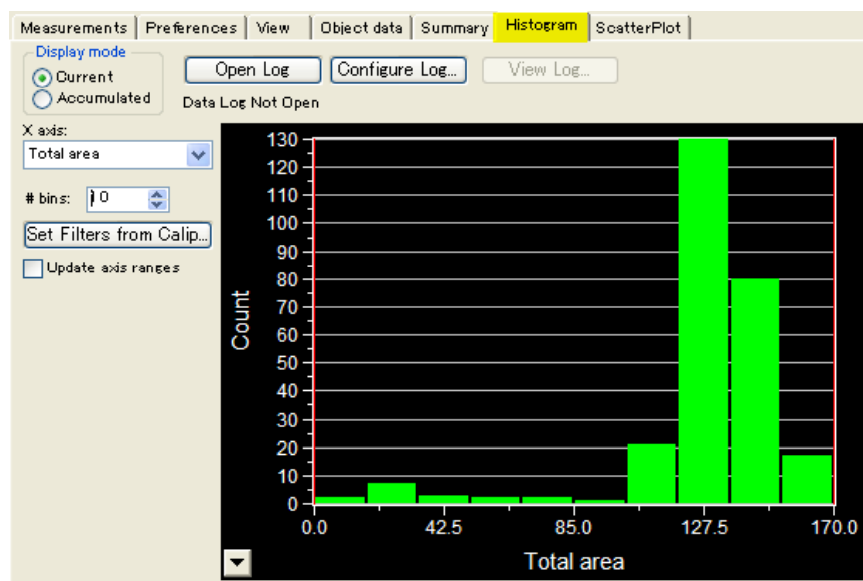
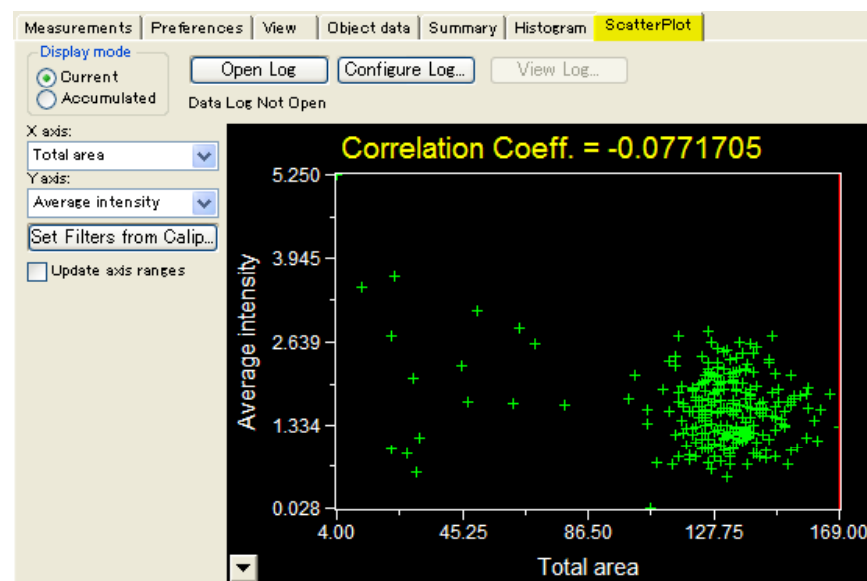
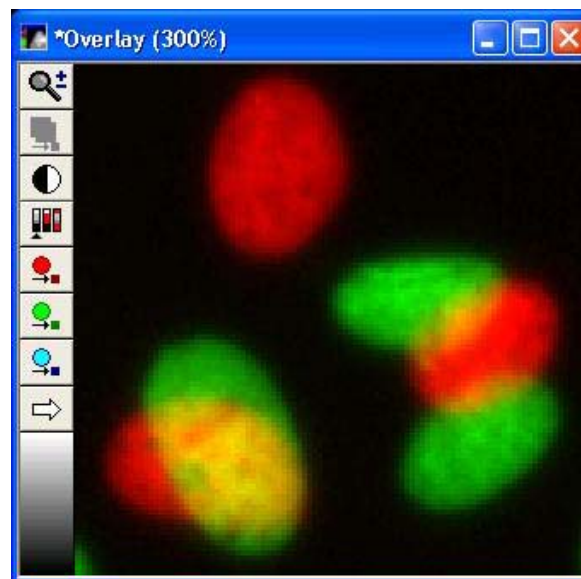
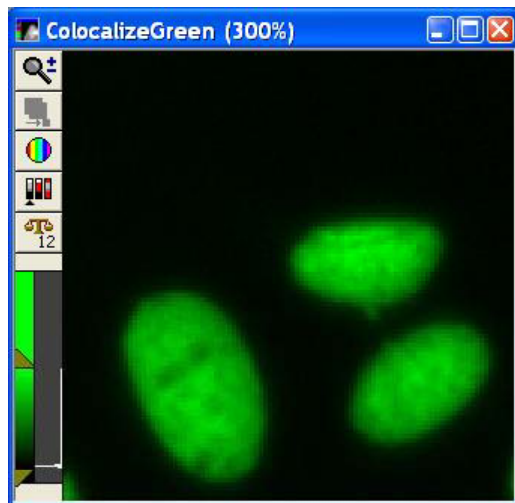
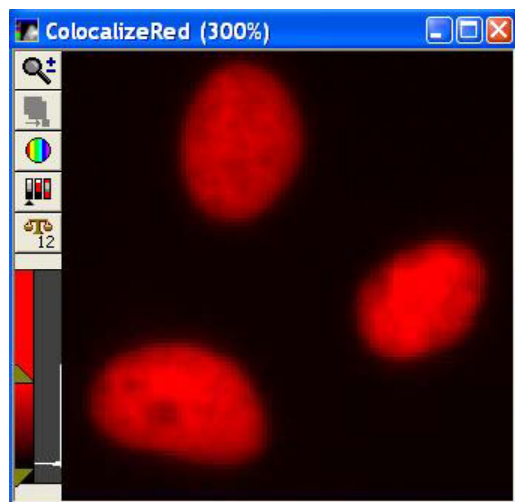
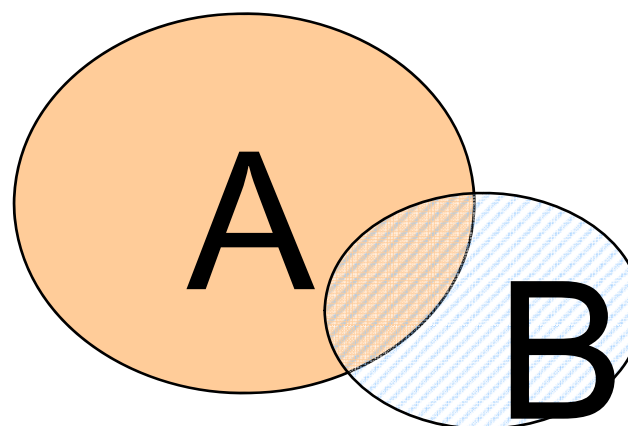
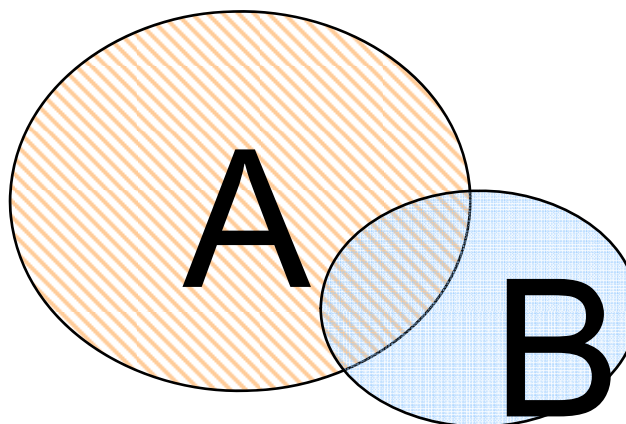
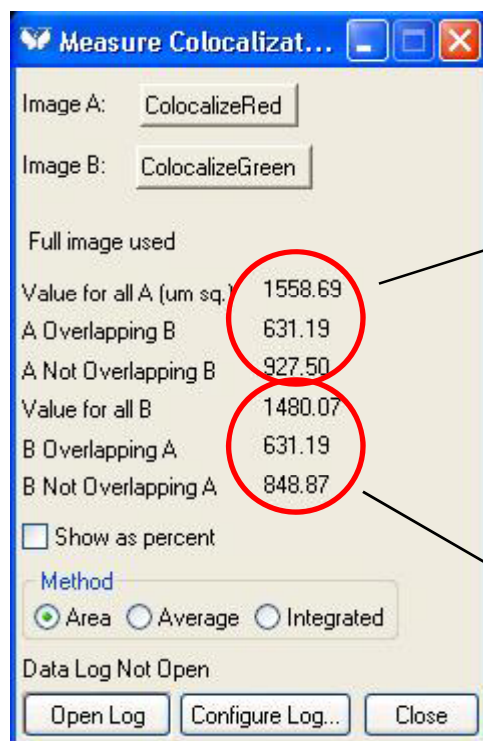
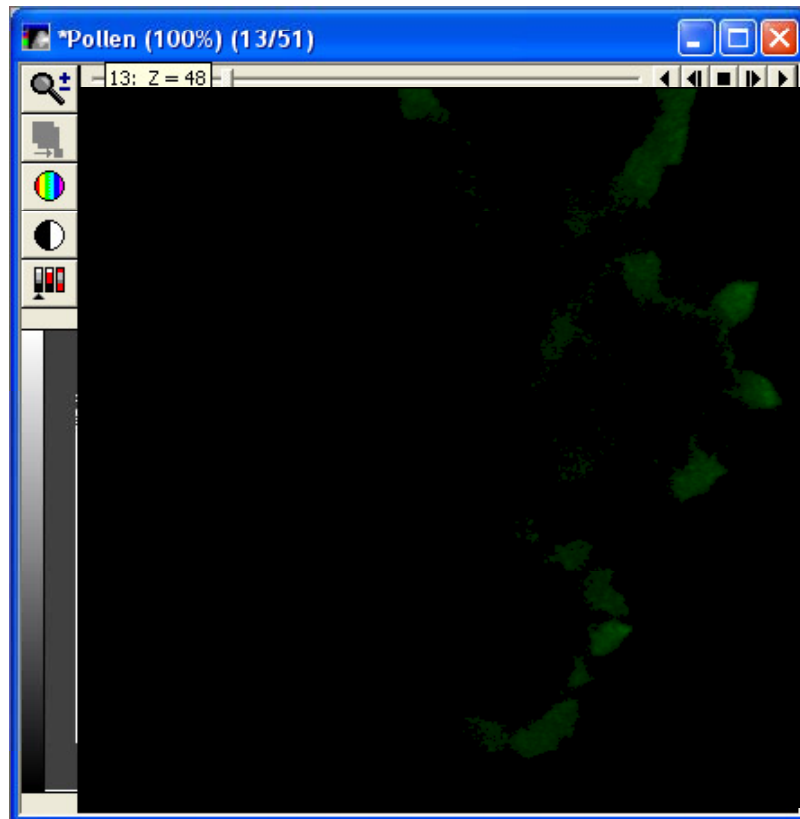
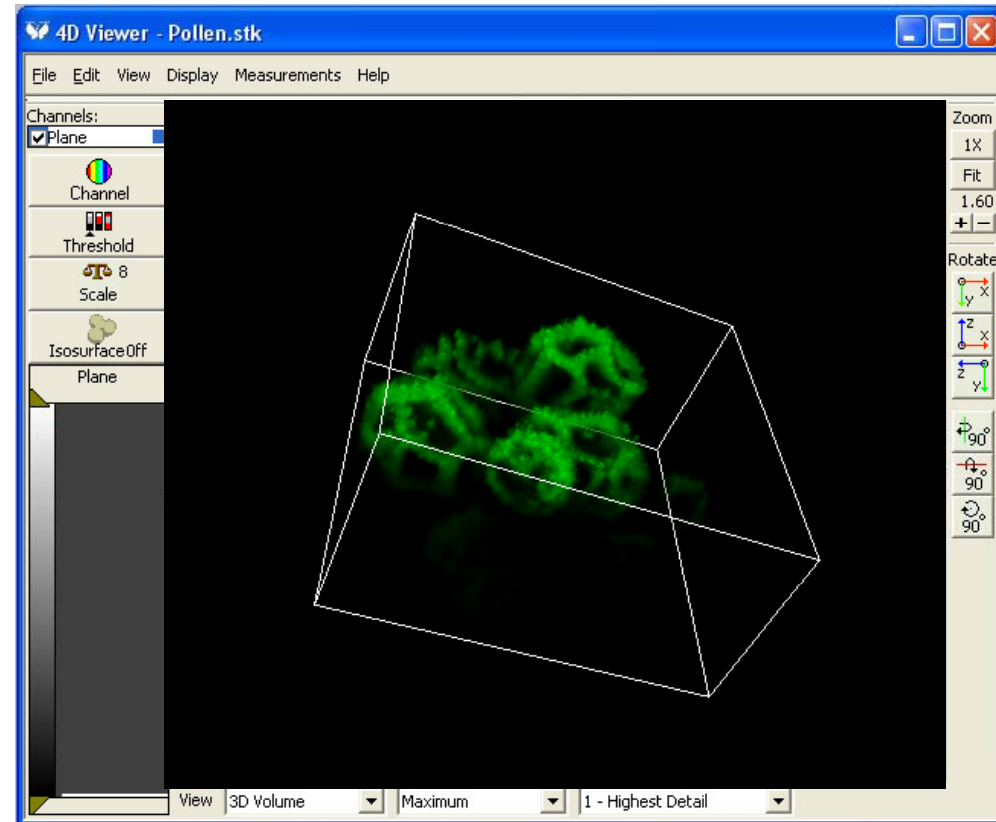
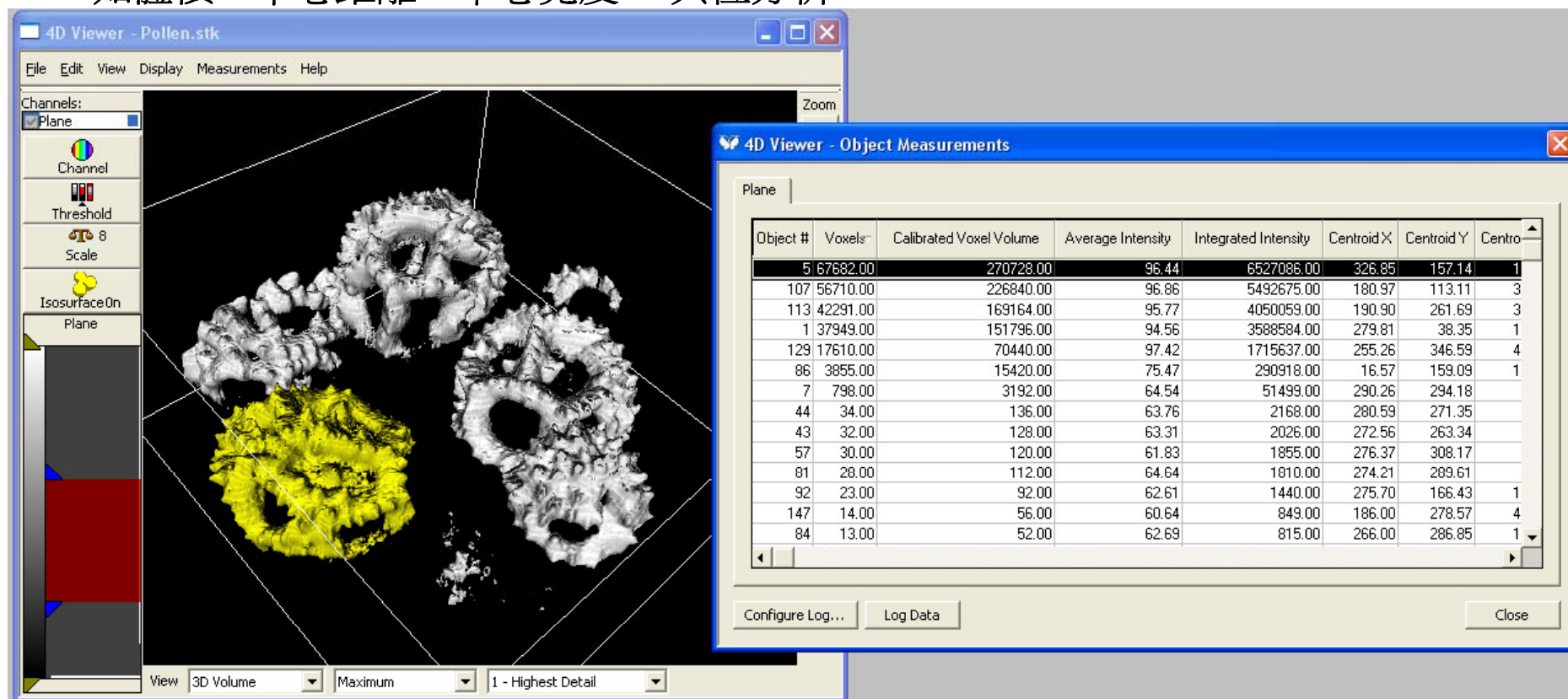
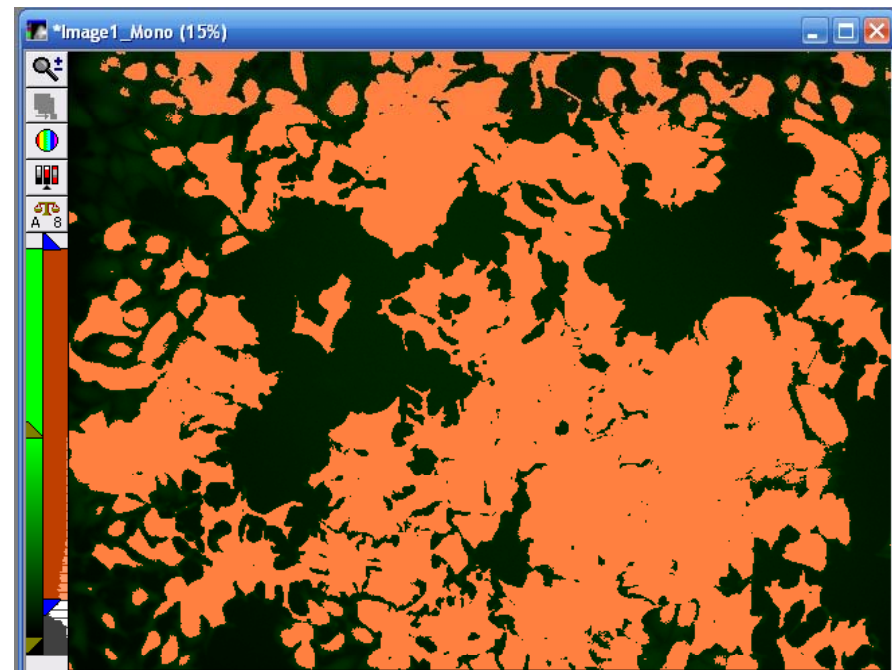
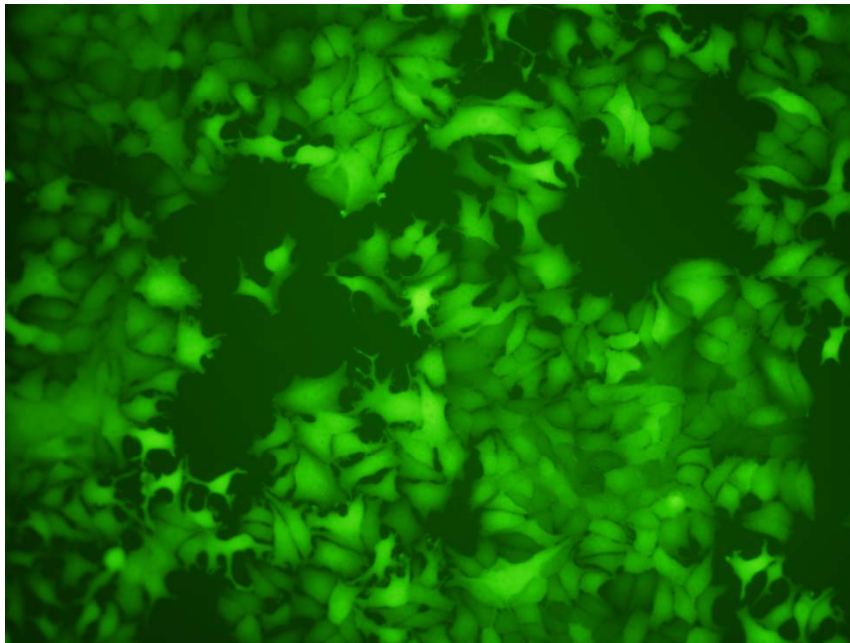
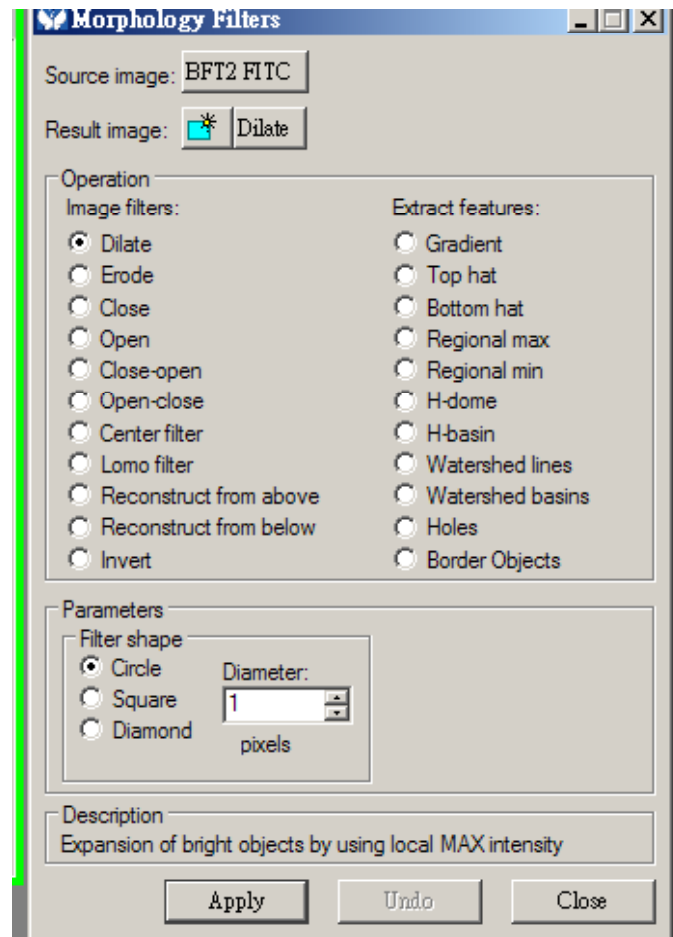
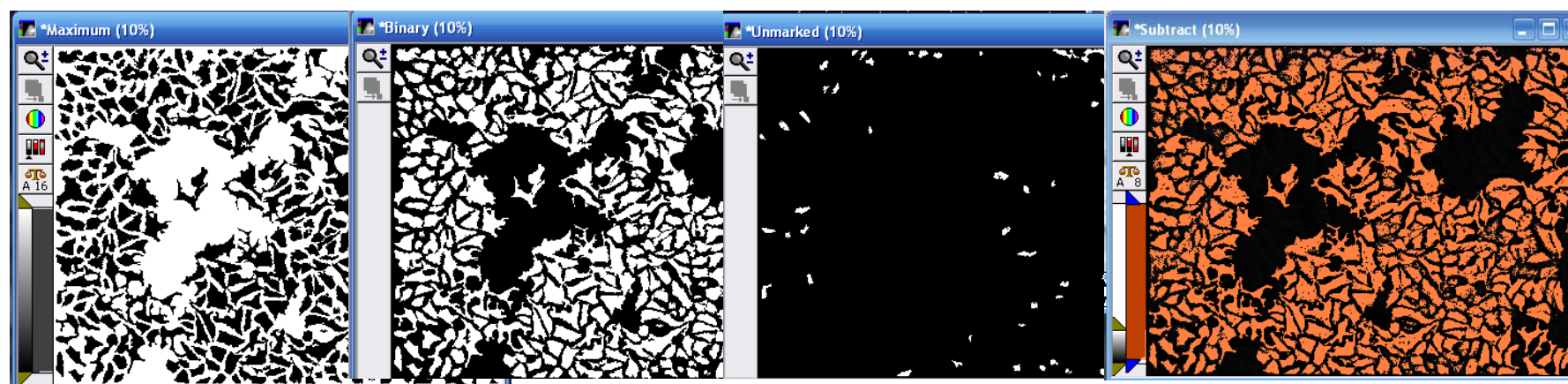
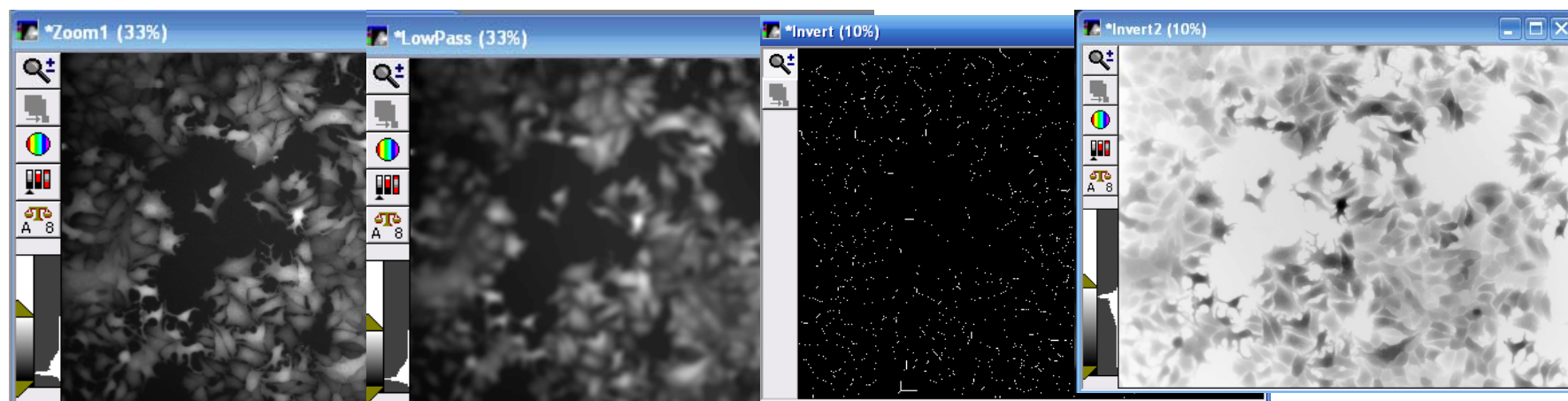
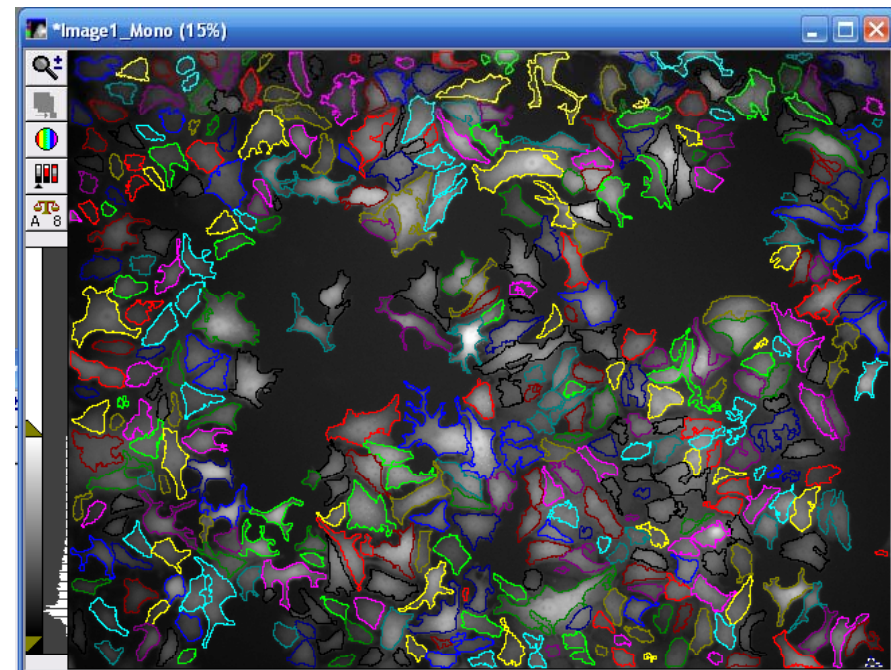
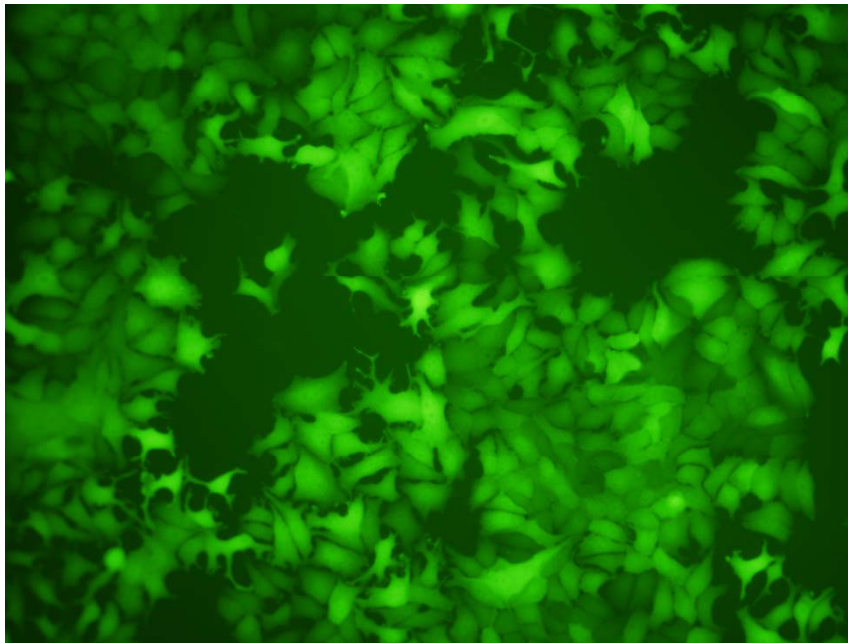
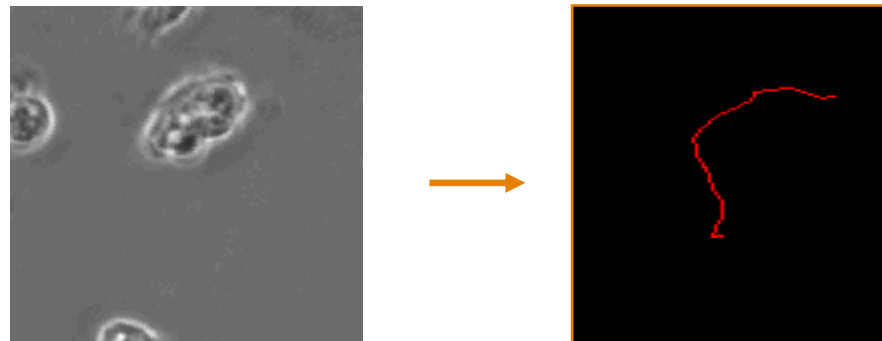
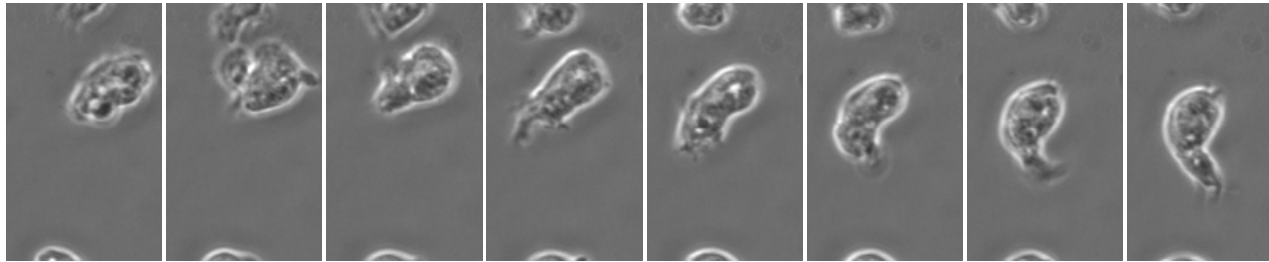
Print Table

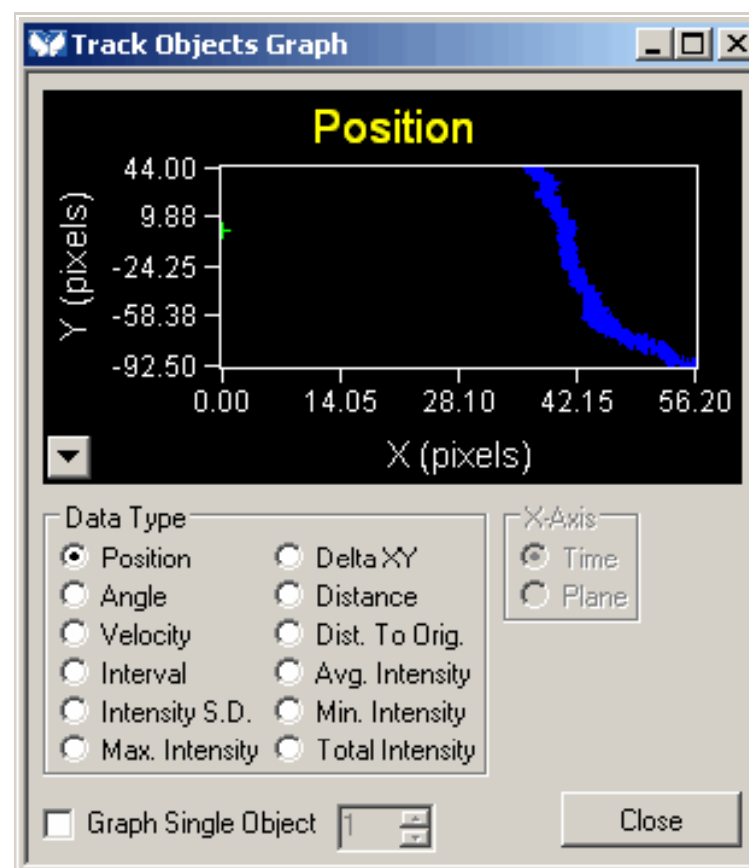Close


Multi Dimensional Motion Analysis

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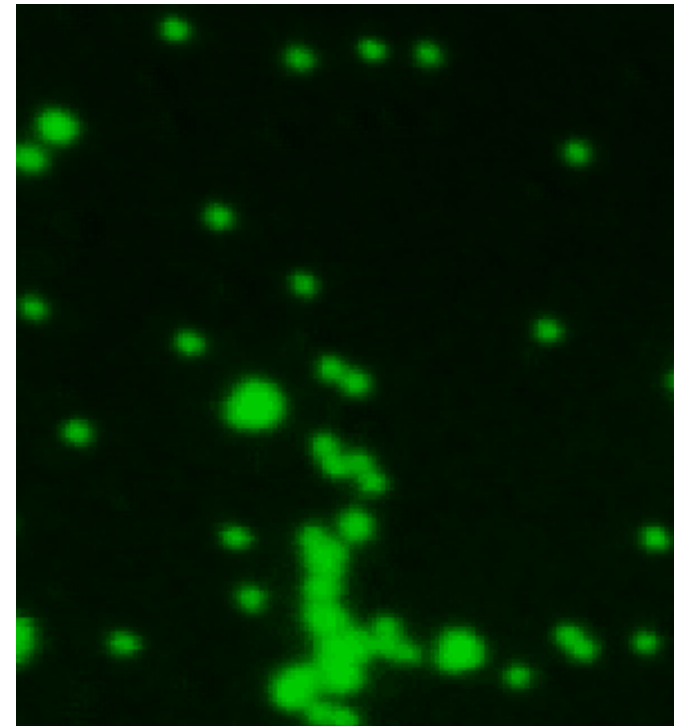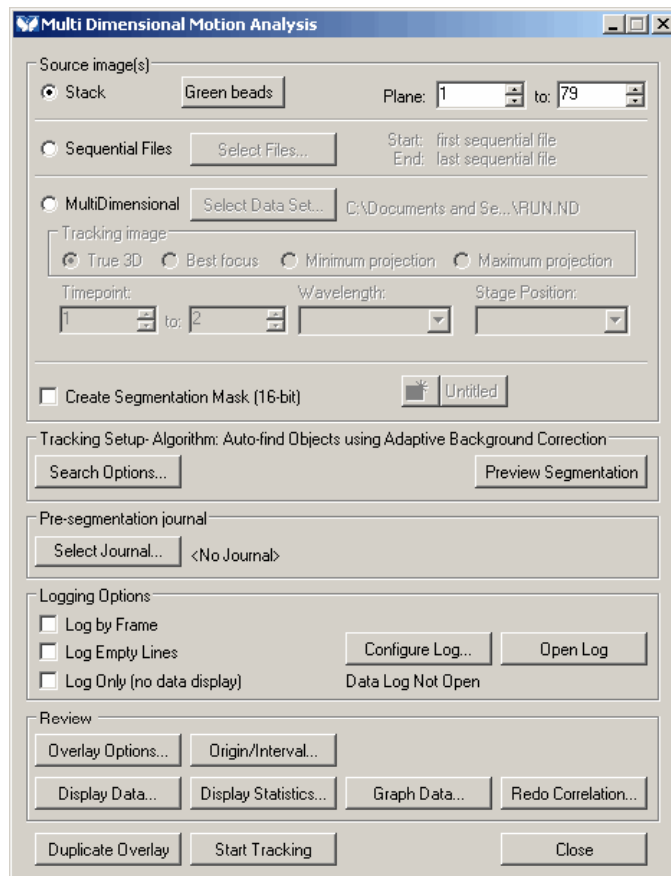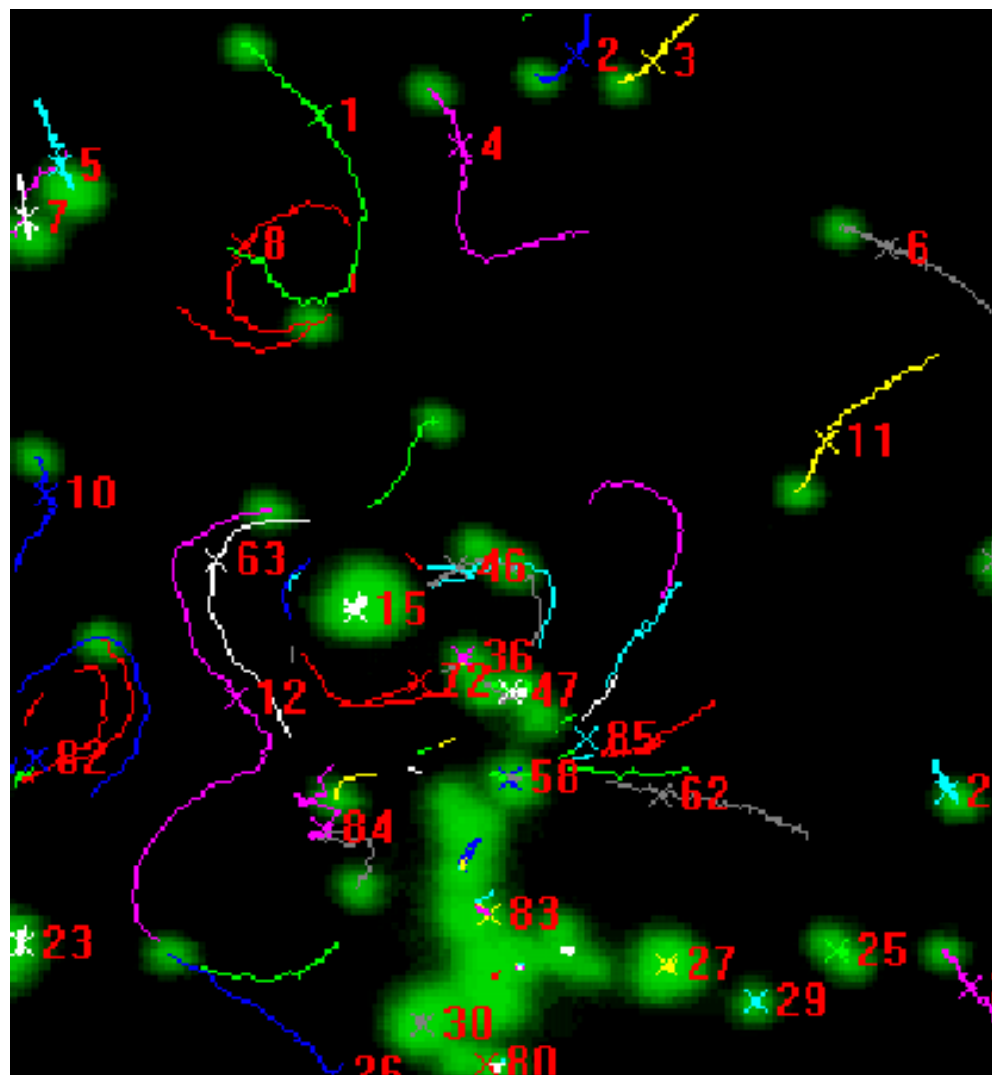
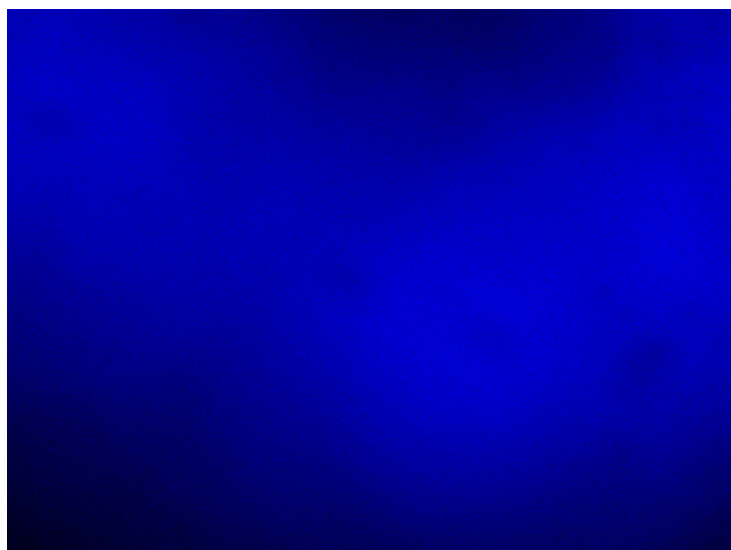
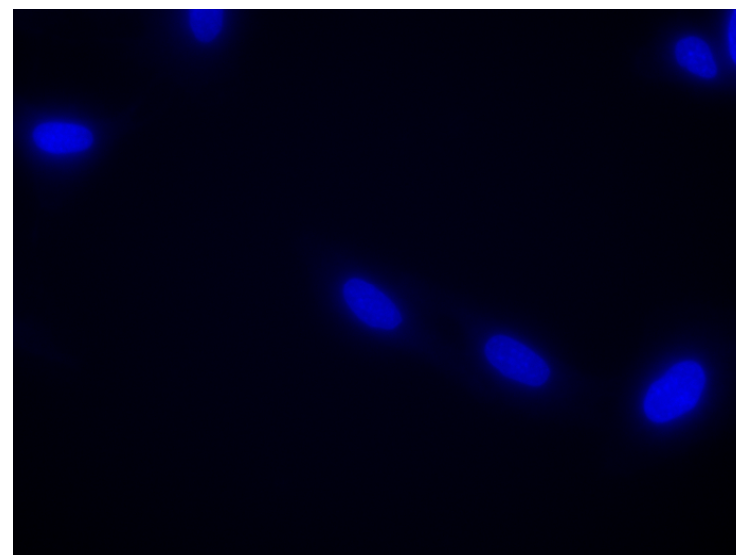
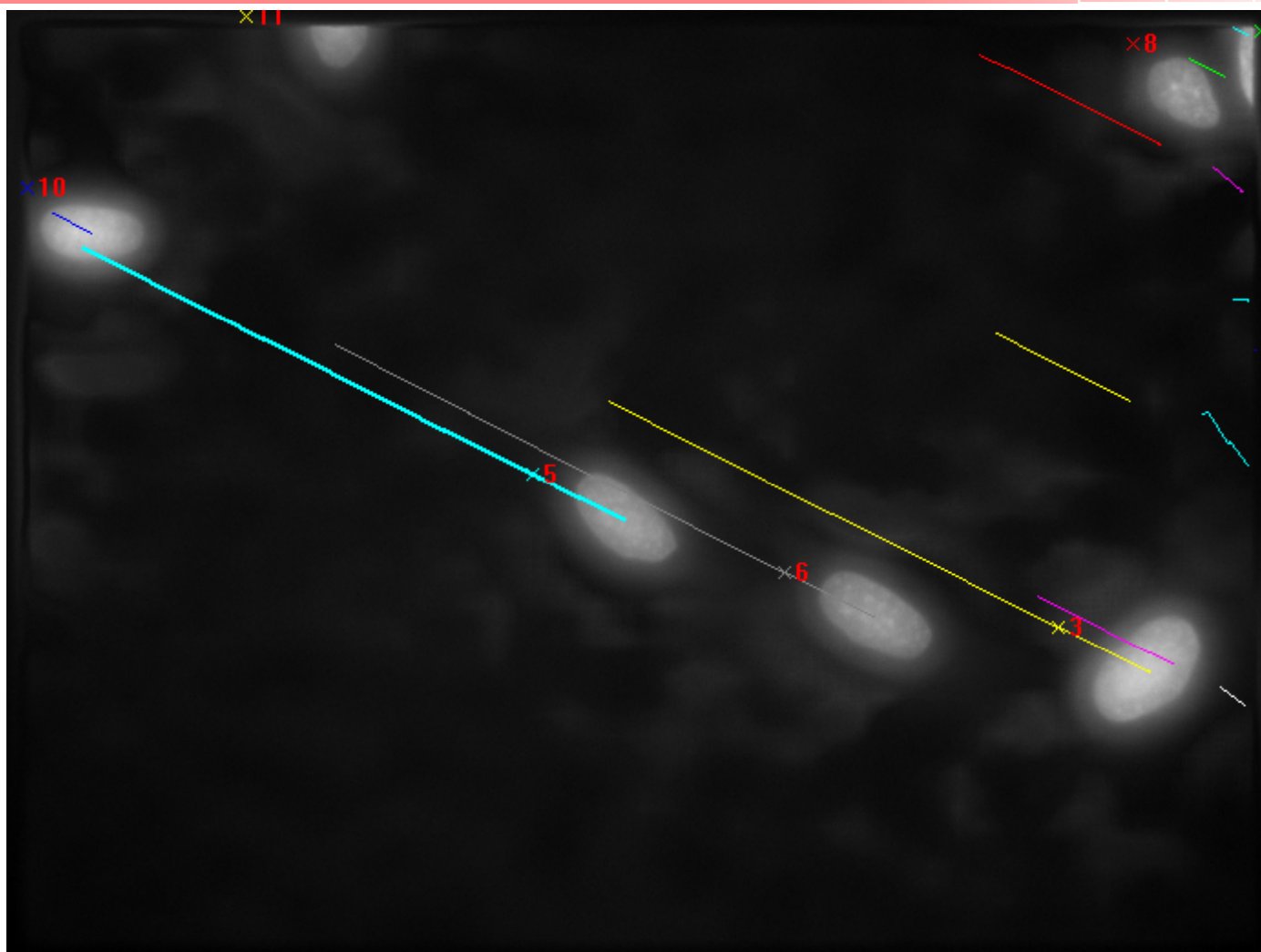
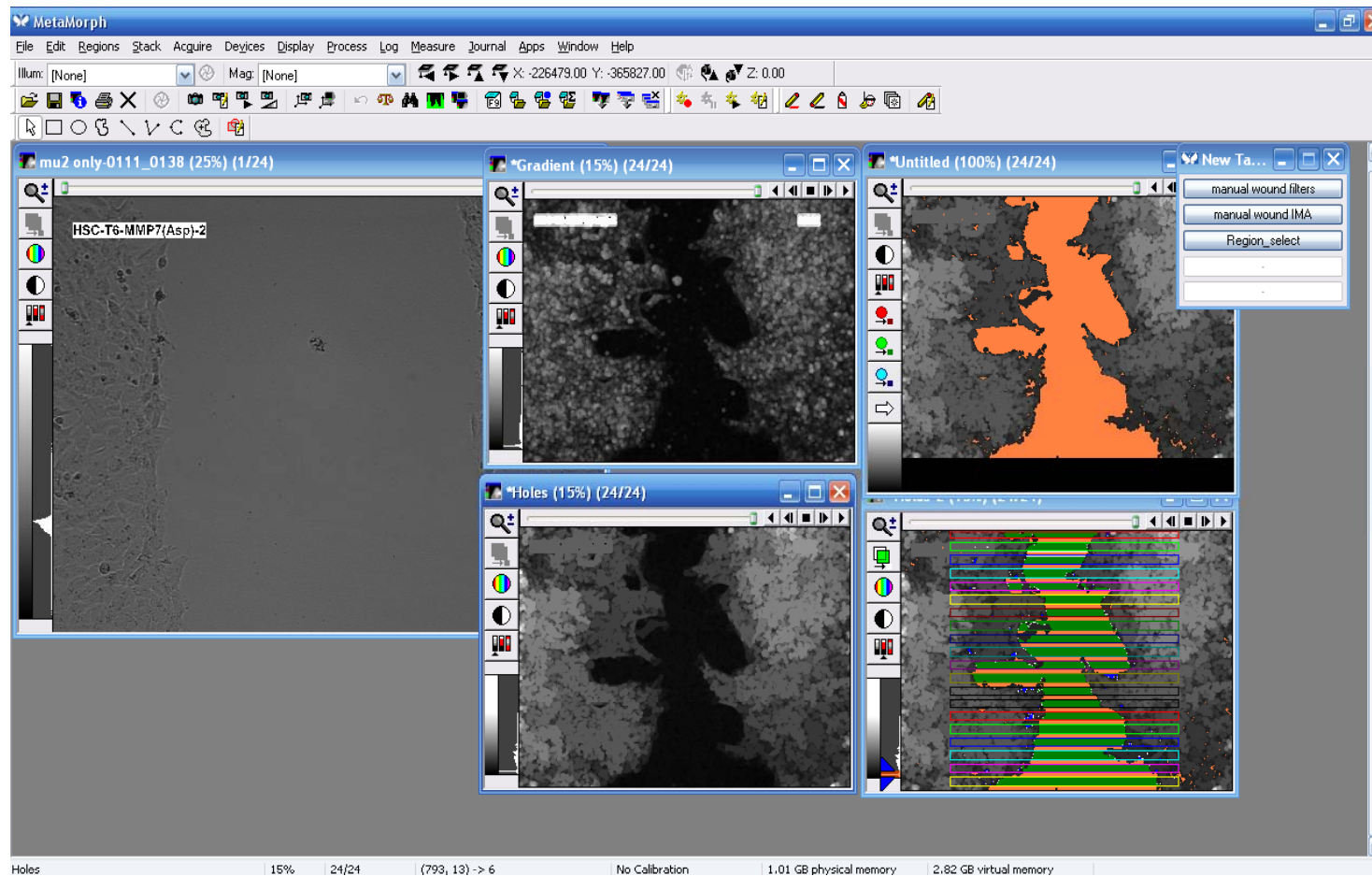
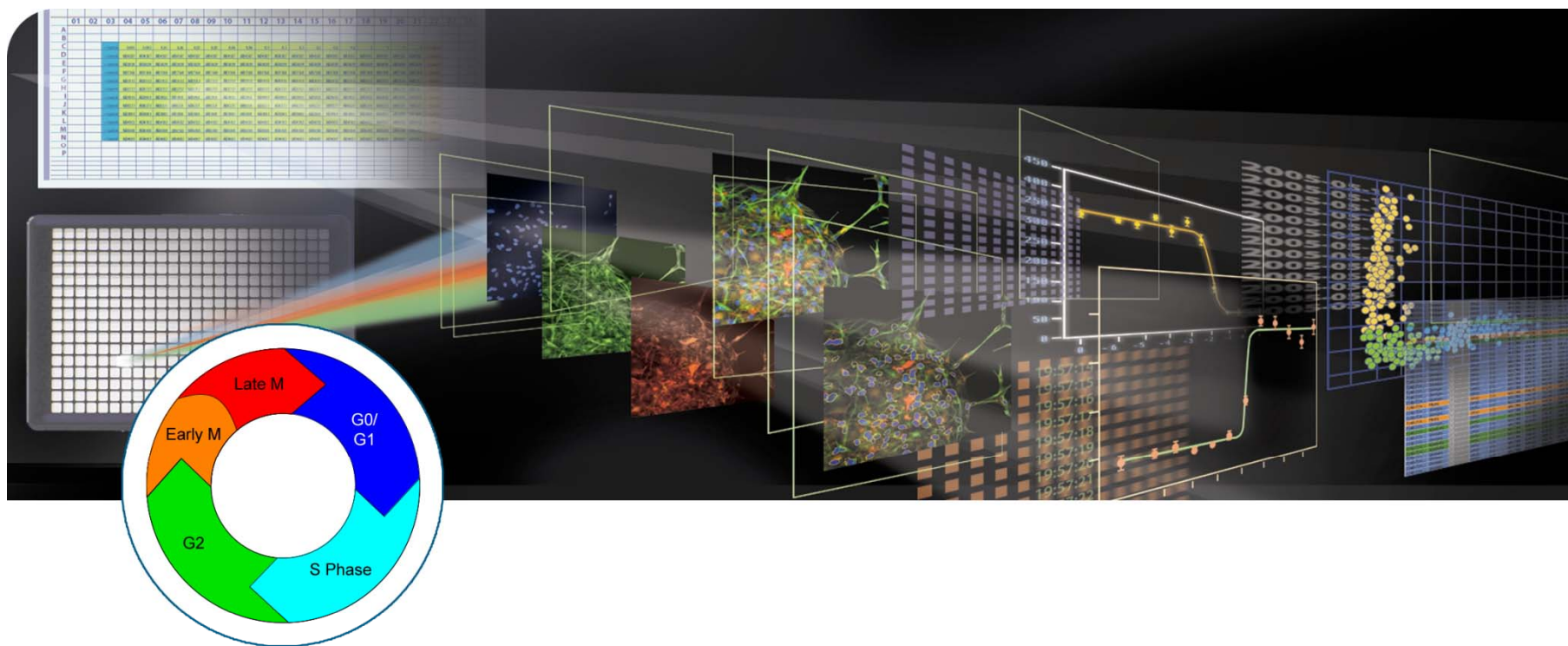
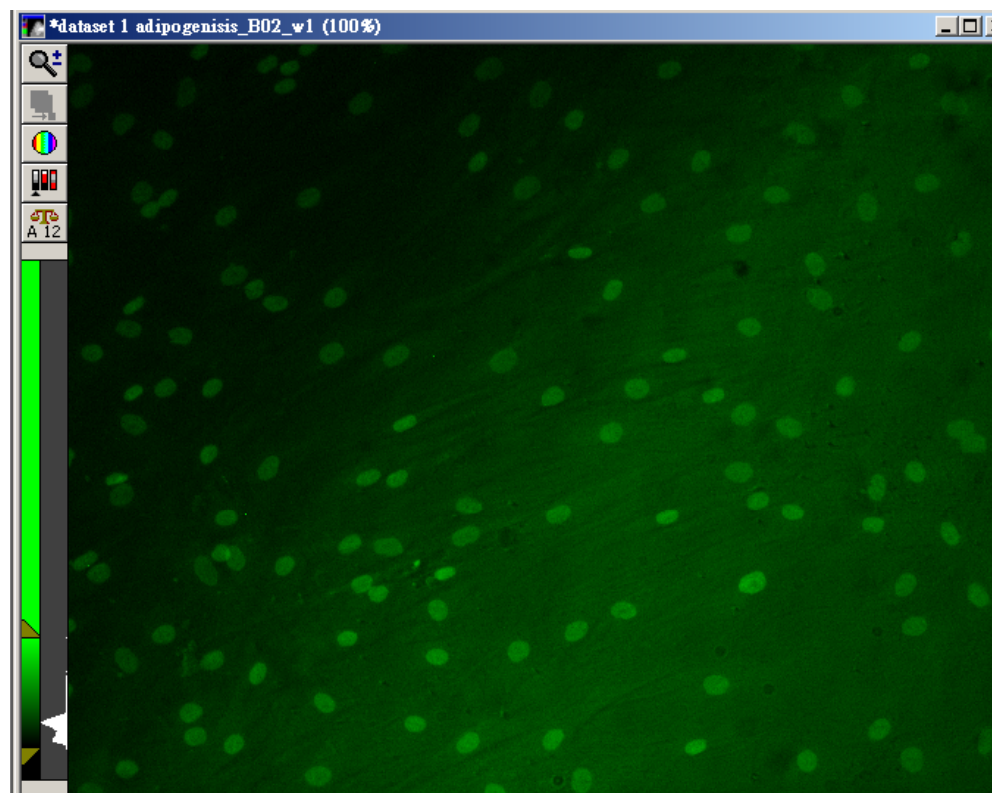
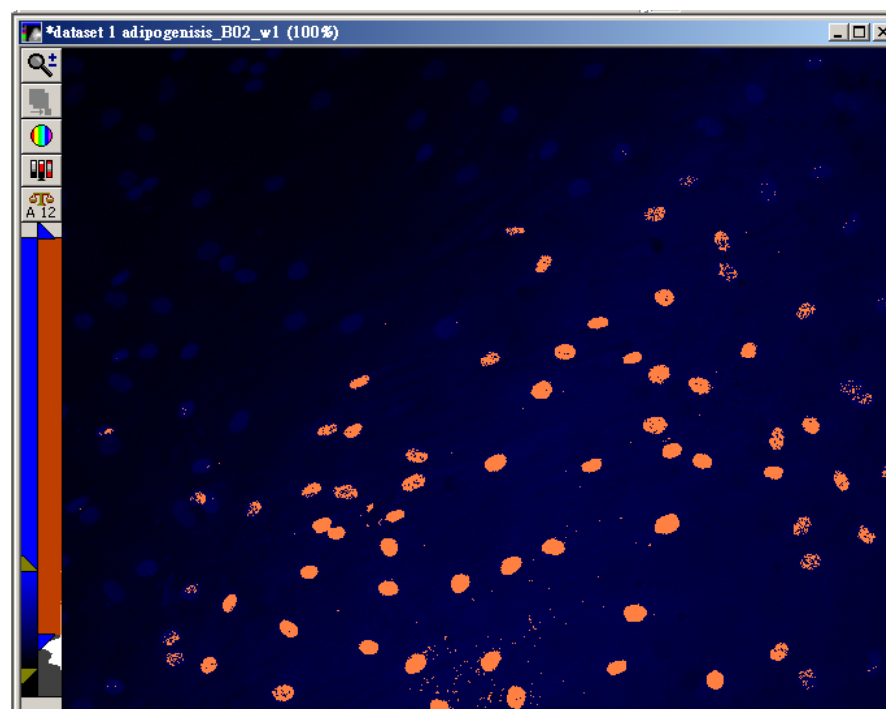
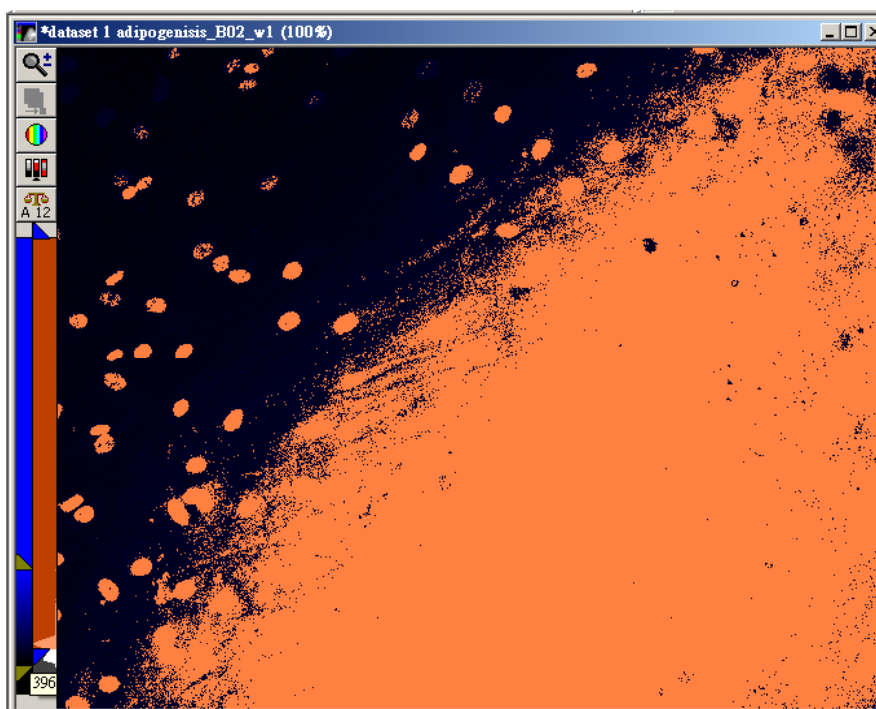
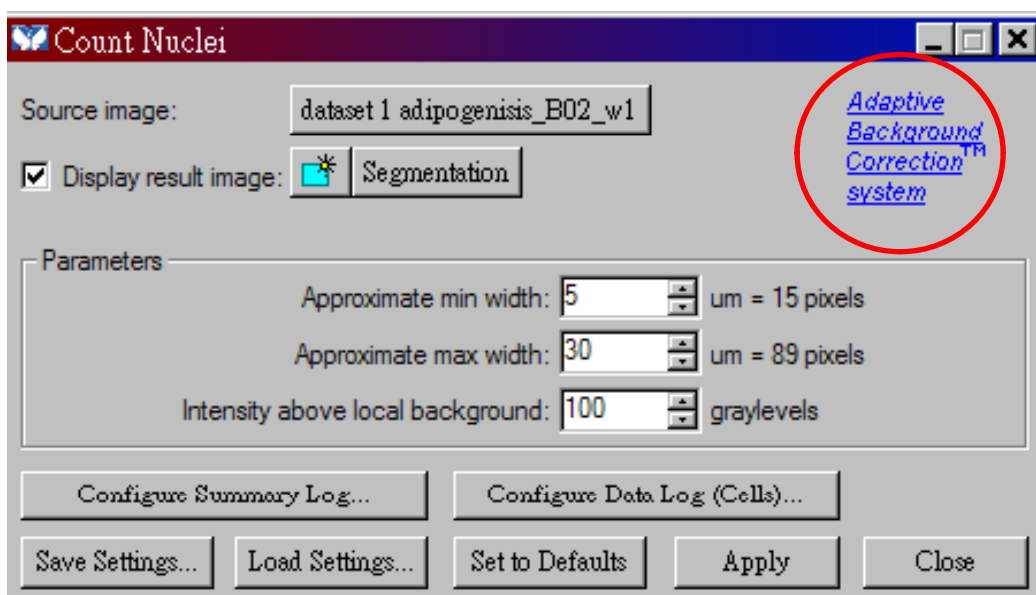
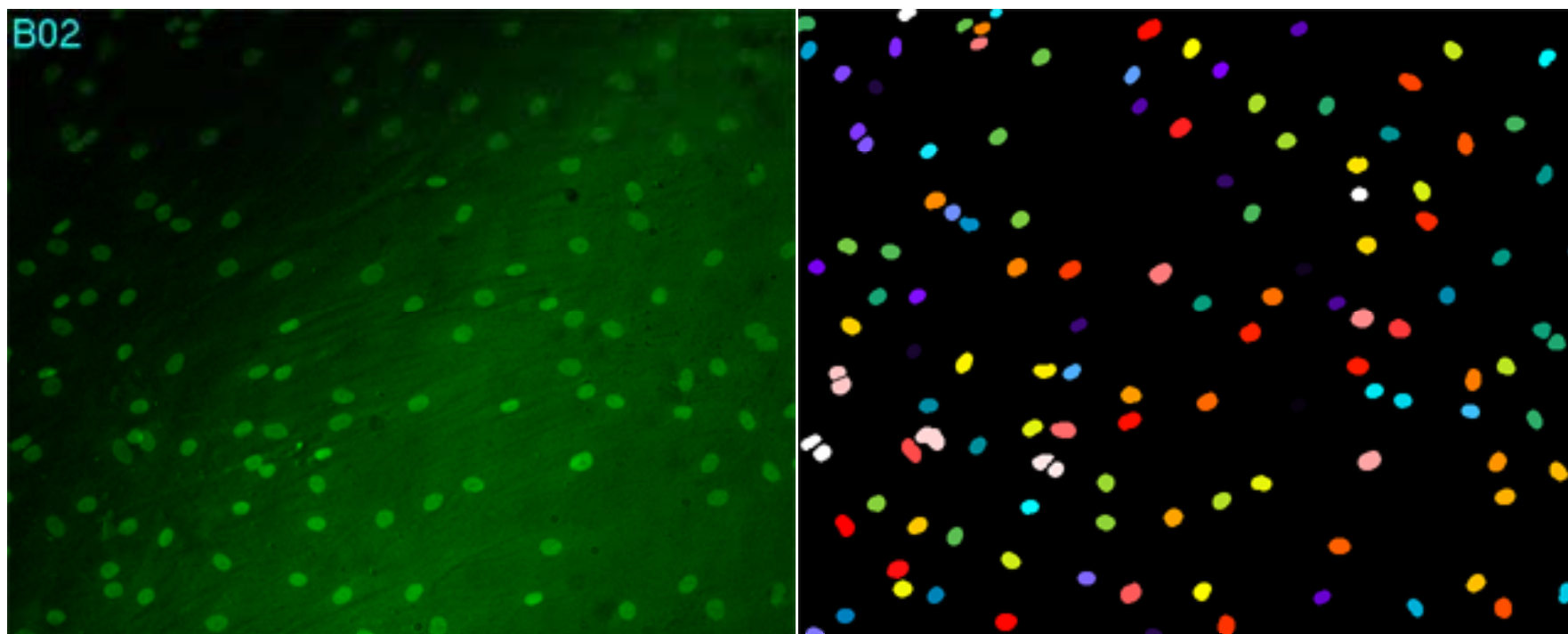
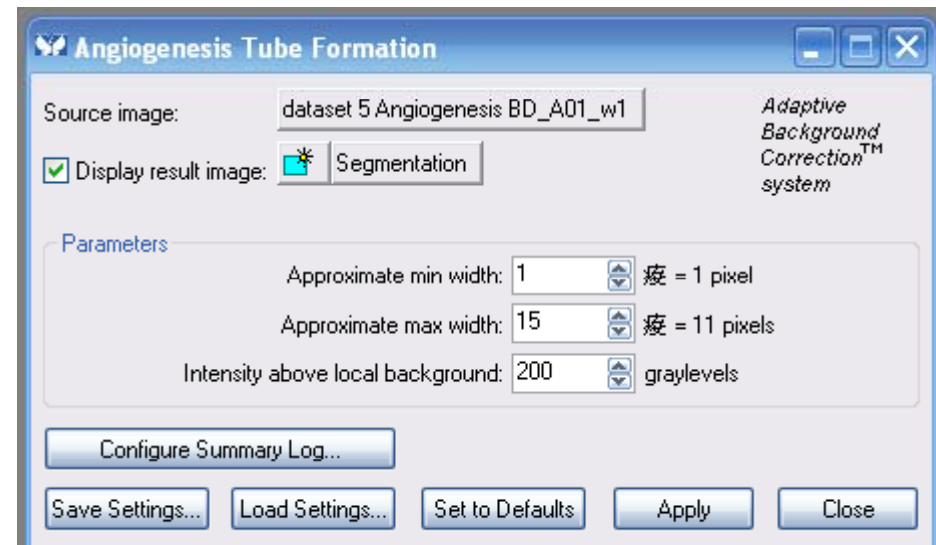
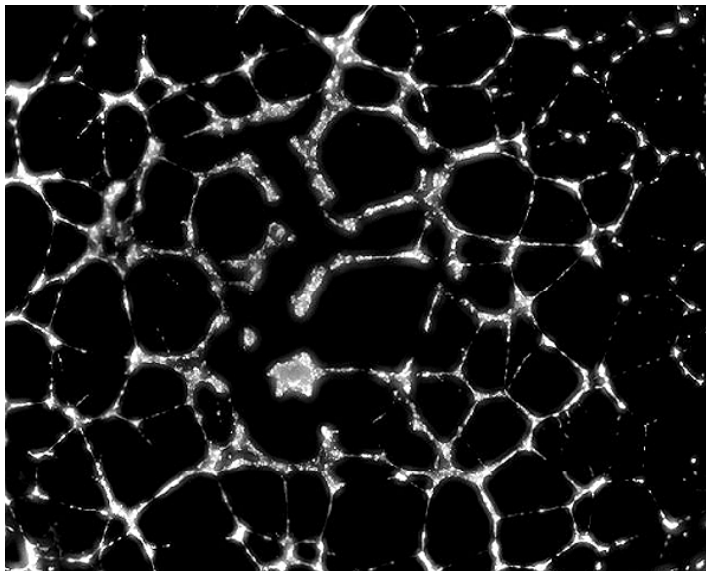
Angiogenesis Results

Angiogenesis Results

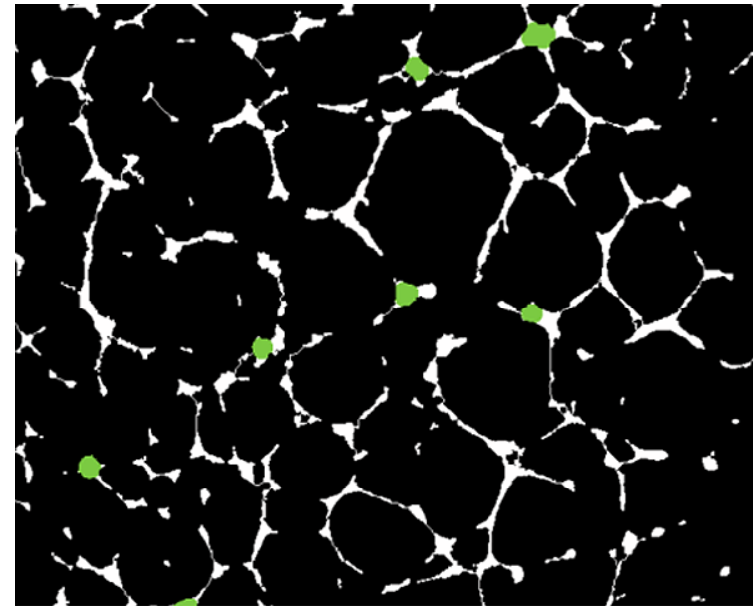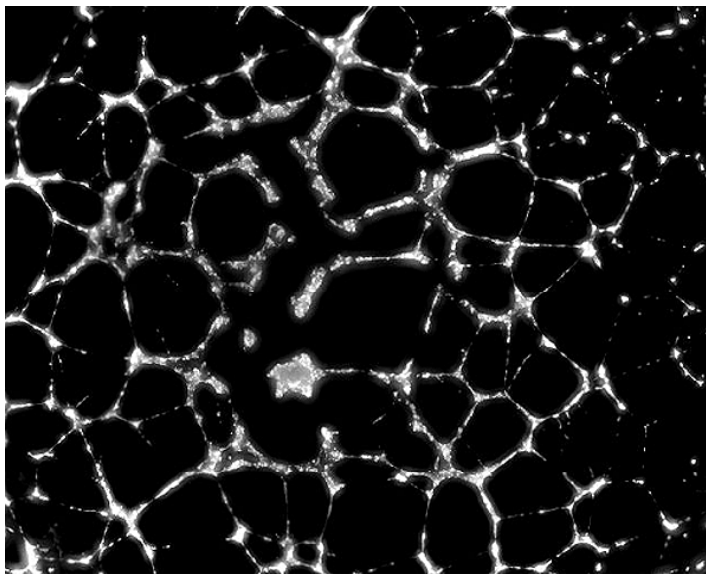Left: from stack, a **best focus** image is obtained. **Right:** analysis. White: tubes, green: nodes. BD BioCoat™ Angiogenesis System: Angiogenesis Endothelial Cell Tube Formation Assay. Courtesy of Min Wu, BD Biosciences.

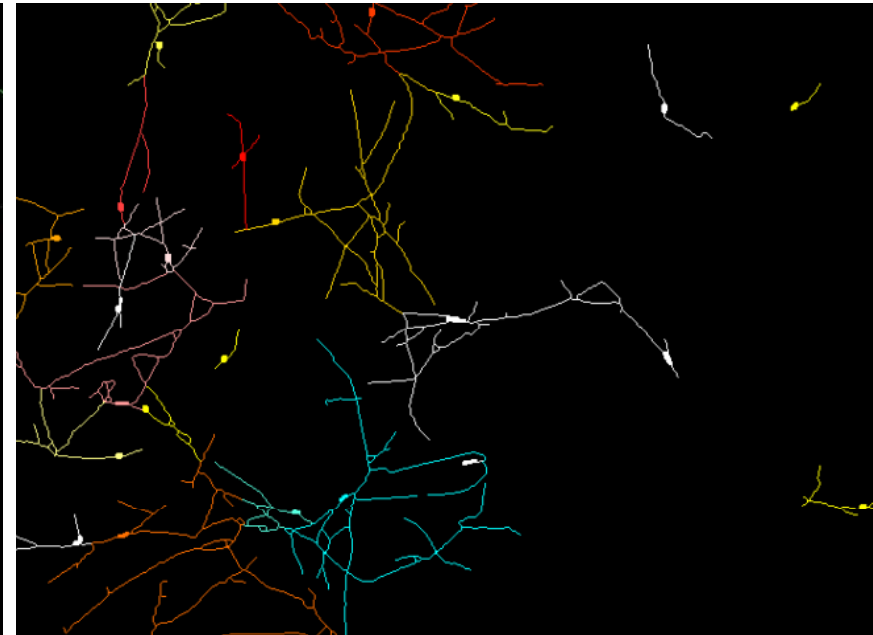
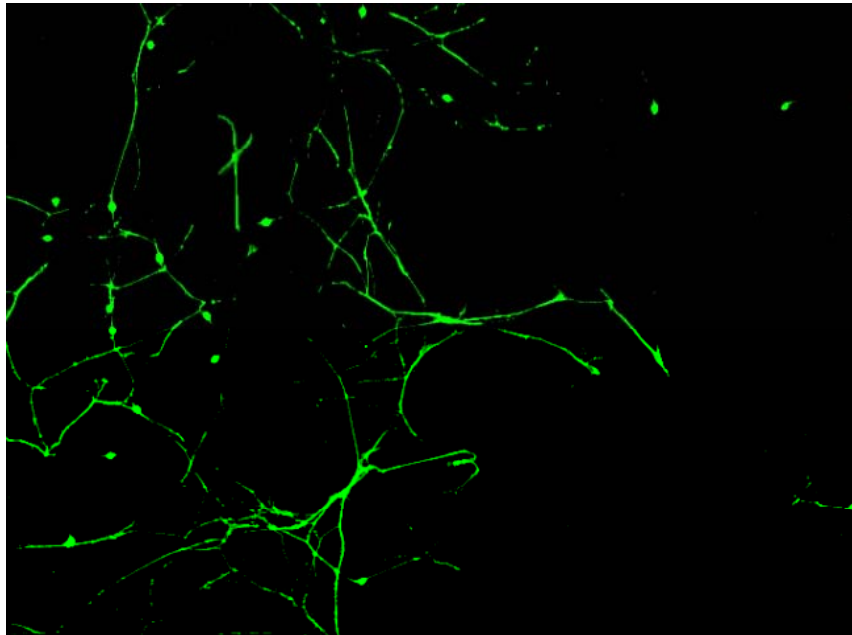
Neurite Outgrowth Application Module

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Neurite Outgrowth Results


Left: image acquisition. Right: Each filament is assigned to a cell body. All the filaments and cell bodies are then measured. Images courtesy of Kris Poulsen and Davide Foletti, Rinat Neuroscience Corporation.

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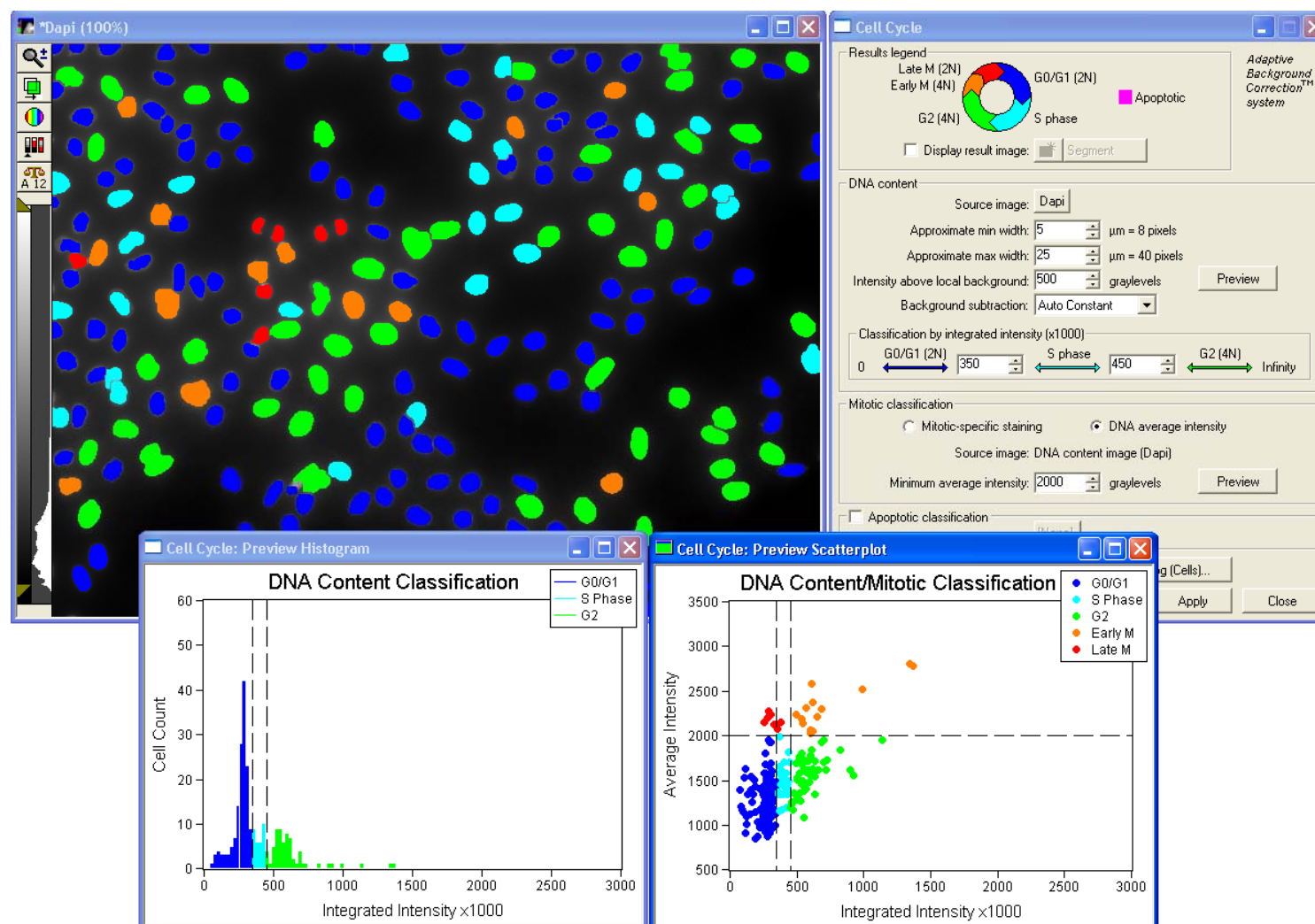
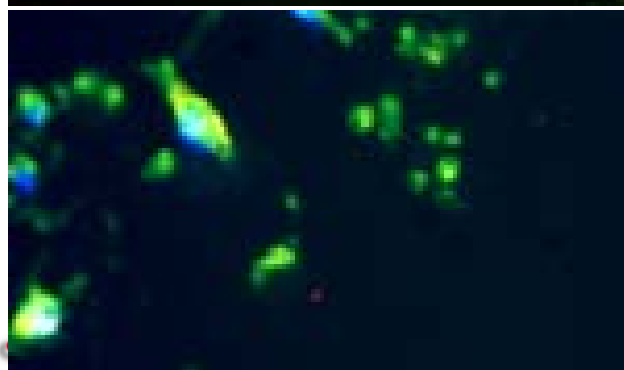
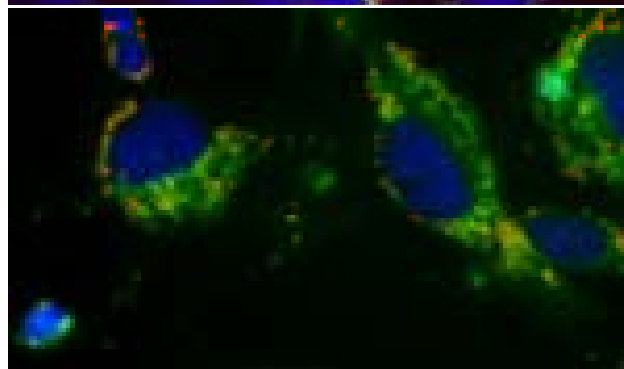
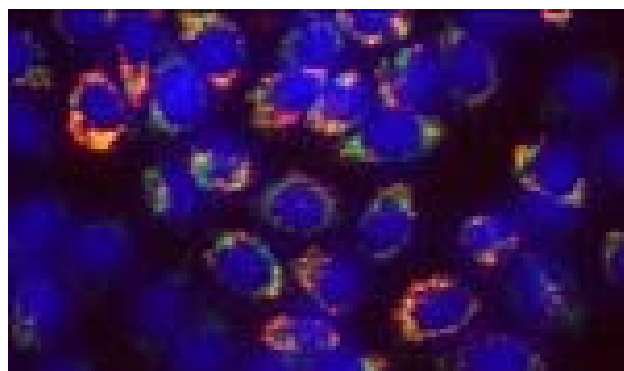
1. Vybrant #7 from Molecular Probes
2. Annexin V
3. JC-1 with Hoechst: used to study loss of mitochondrial potential

Cell Health Results

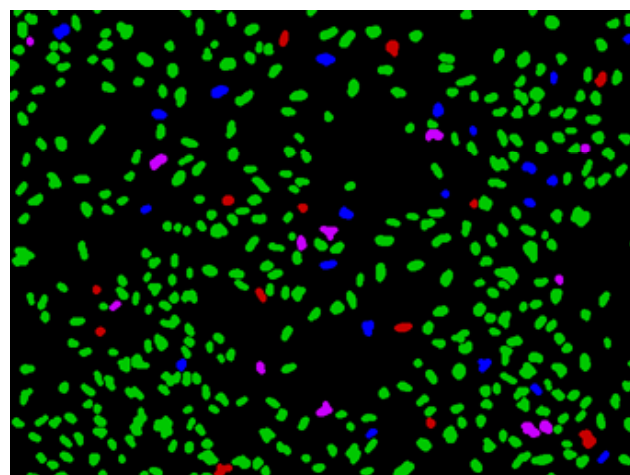Adherent Chinese Hamster Ovary (CHO-K1) cells were incubated with various concentrations of staurosporine. From top to bottom: control, 0.1 μM staurosporine, and 3 μM staurosporine. BIOMOL International's Mit-E-Y Mitochondrial Permeability Detection (Hoechst 33342 + JC-1). Courtesy of BIOMOL International.

Cells analyzed by the module. Green: viable, blue: early apoptotic, purple: late apoptotic, red: necrotic.

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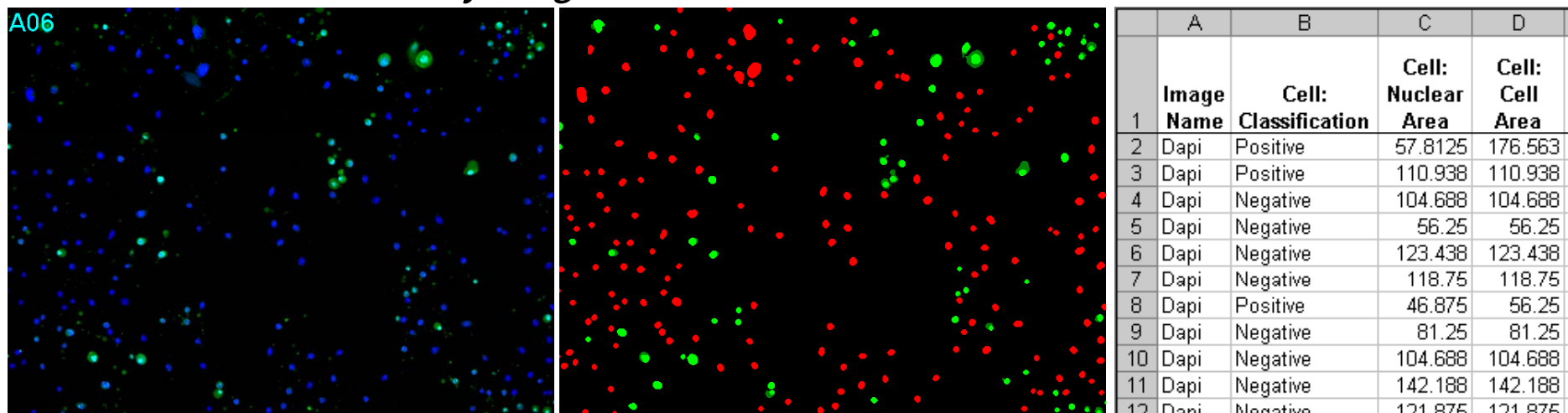
Left: Acquired image, two wavelengths overlay. **Center:** Analyzed image shows the identification of two probes. **Right:** Data collected can be exported to a spreadsheet program for further analysis.

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Major Instruments Co., Ltd.